English

Installation instructions

Weighing Table XP26PC/XPE26PC

Deutsch

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
Wägetisch XP26PC/XPE26PC

Français

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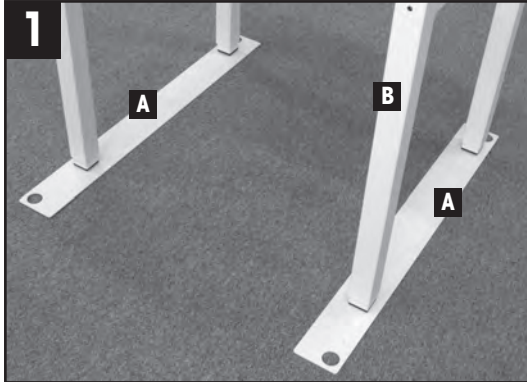
Table de pesage XP26PC/XPE26PC

Choose an appropriate location (no vibrations, see balance manual). Locate positioning guides [A] and place balance table [B] (inner table) into the respective holes.

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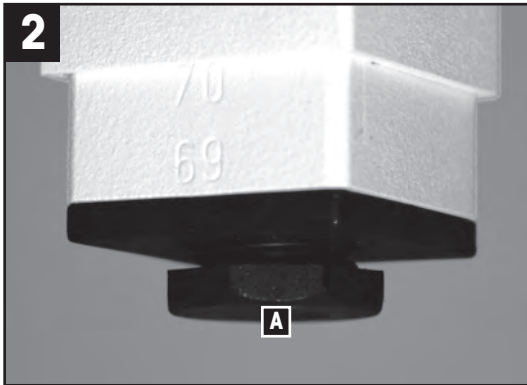
Use the leveling feet [A] to adjust the height of the balance table until the positioning guides can be moved freely (approx. 3mm).

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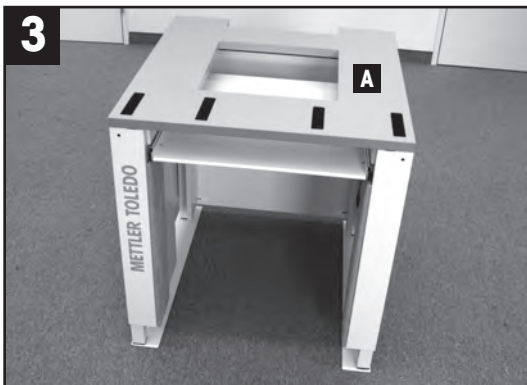Put the working table [A] (outer table) over the balance table, onto the positioning guides.

Caution: This step requires at least two persons!

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
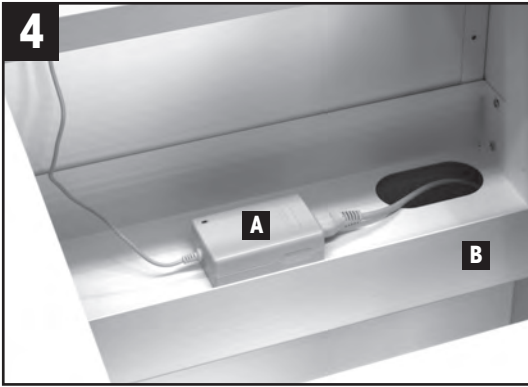
METTLER TOLEDO

Place the AC adapter [A] in the cable tray [B] at the back of the working table.

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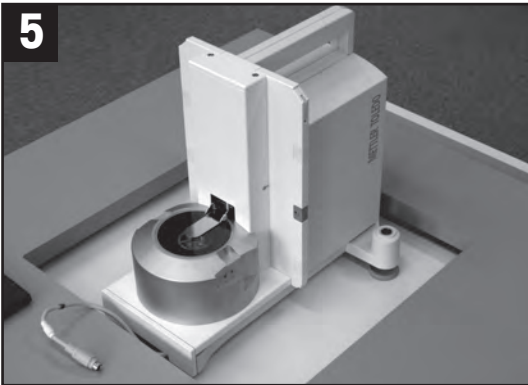
Remove terminal and light barrier from the balance. Place the balance onto the balance table and level it.

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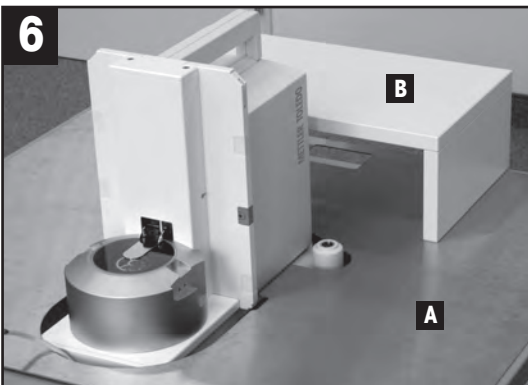
Carefully place the metal sheet [A], then install the console [B]: Insert the two pins at the bottom of the console into the respective slots of the metal sheet (the console fixes the metal sheet).

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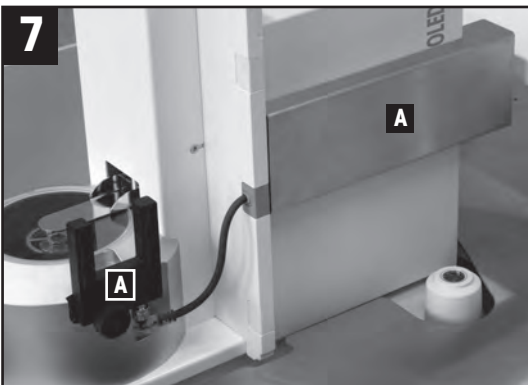
Install the light barrier unit [A] and connect the terminal. Connect the balance to the AC adapter.

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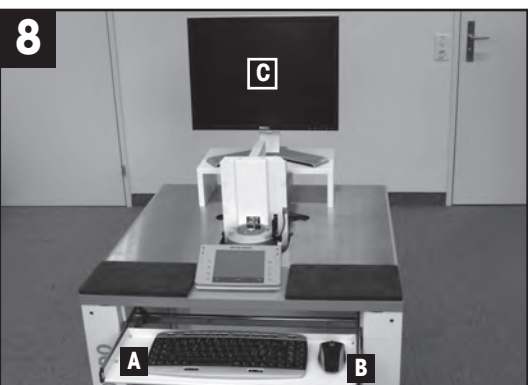
Install wireless keyboard [A] and wireless mouse [B]. Install display [C], if available.

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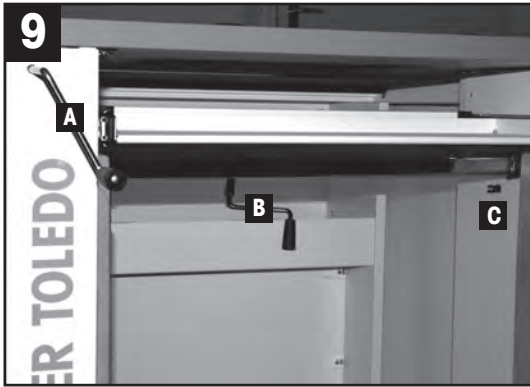
Use the two handles [A and B] to adjust both tables to an optimum ergonomic height. Then, clip the handle [A] into the respective holder [C].

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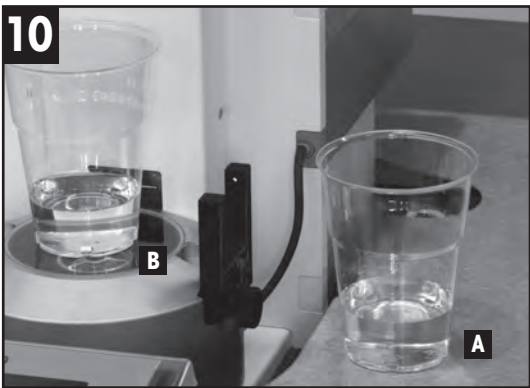
Check installation with two cups of water: Slightly shake the working table [A]: The water surface on the balance table [B] should remain still.

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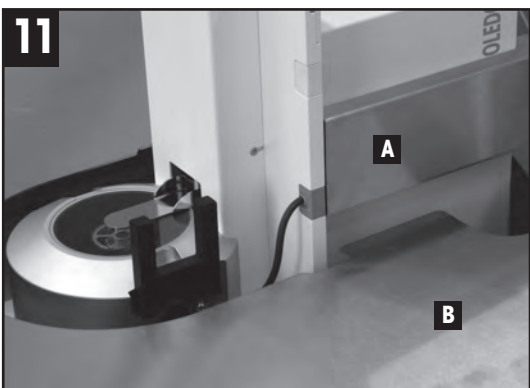CAUTION: When adjusting the height of the tables, the cover [A] and the metal sheet [B] must not touch, otherwise the balance might get damaged.

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© Mettler-Toledo AG 12 /2014
Mettler-Toledo AG, Laboratory Weighing