English

Installation instructions

Weighing Table XP26/XPE26, XP56/XPE56

Deutsch

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
Wägetisch XP26/XPE26, XP56/XPE56

Français

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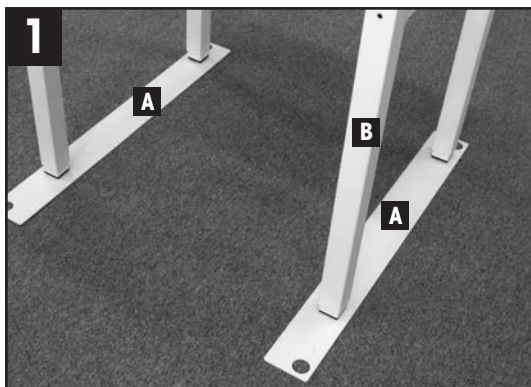
Table de pesage XP26/XPE26, XP56/XPE56

Choose an appropriate location (no vibrations, see balance manual). Locate positioning guides [A] and place balance table [B] (inner table) into the respective holes.

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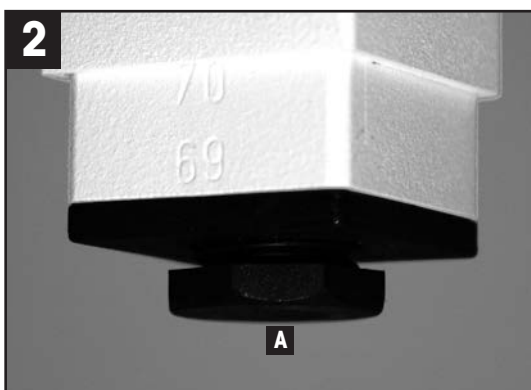Use the leveling feet [A] to adjust the height of the balance table until the positioning guides can be moved freely (approx. 3mm).

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Put the working table [A] (outer table) over the balance table, onto the positioning guides.

Caution: This step requires at least two persons.

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
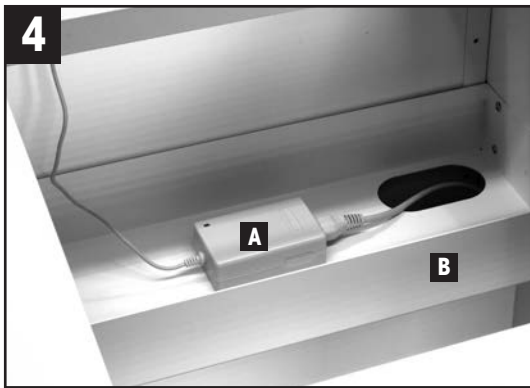
METTLER TOLEDO

Place the AC adapter [A] in the cable tray [B] at the back of the working table.

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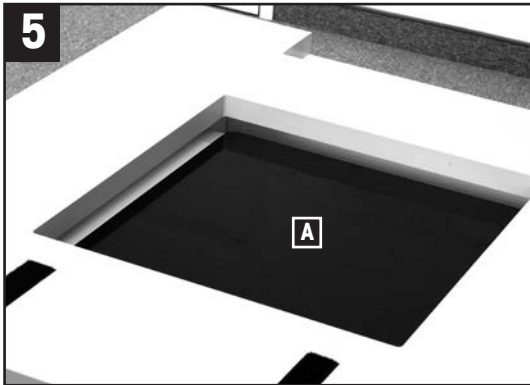
Carefully place the stone plate [A] in the balance table.

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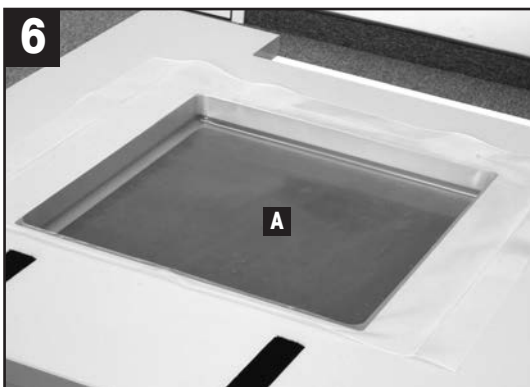
Place the flexible drip pan [A] in the working table. Adjust the height of the balance table (see step 5) so that the bottom of the drip pan is properly seated on the stone plate.

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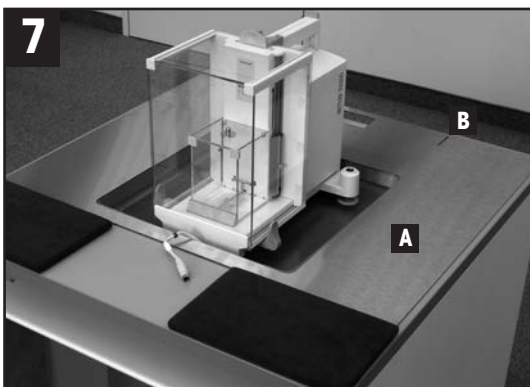
Place the balance, level it and connect it to the AC adapter. Place the metal sheet [A] and align the slots [B] with the recess in the working table.

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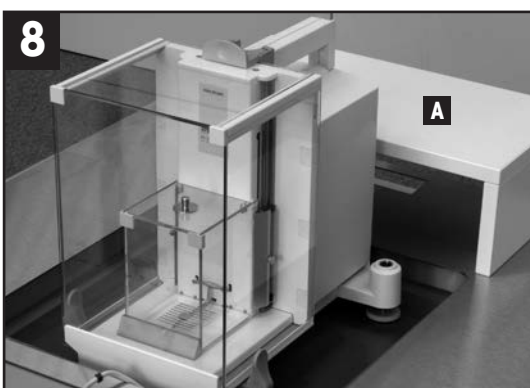
Install console [A]: Insert the two pins at the bottom of the console into the respective slots of the metal sheet (the console secures the metal sheet).

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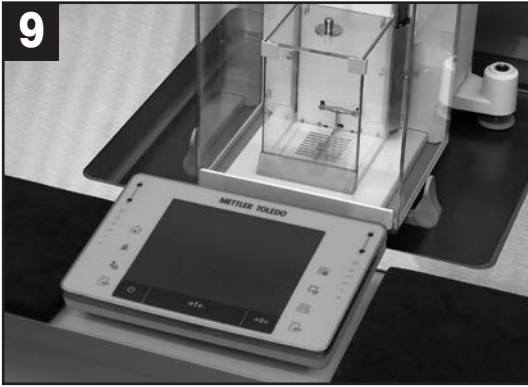Terminal anschliessen.

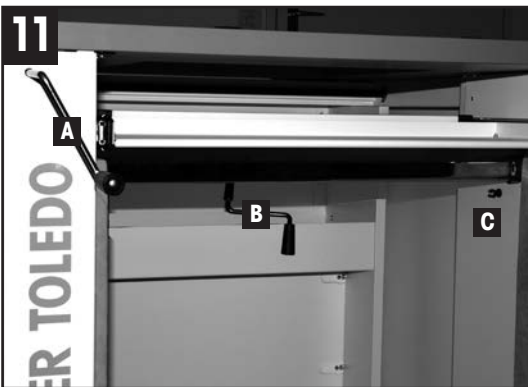
Brancher le terminal.

Install peripheral devices if available.

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Use the two handles [A and B] to adjust both tables to an optimum ergonomic height. Then, clip the handle [A] into the respective holder [C].

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Check installation with two cups of water: Slightly shake the working table [A]: The water surface on the balance table [B] should remain still.

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© Mettler-Toledo AG 12/2014
Mettler-Toledo AG, Laboratory Weighing