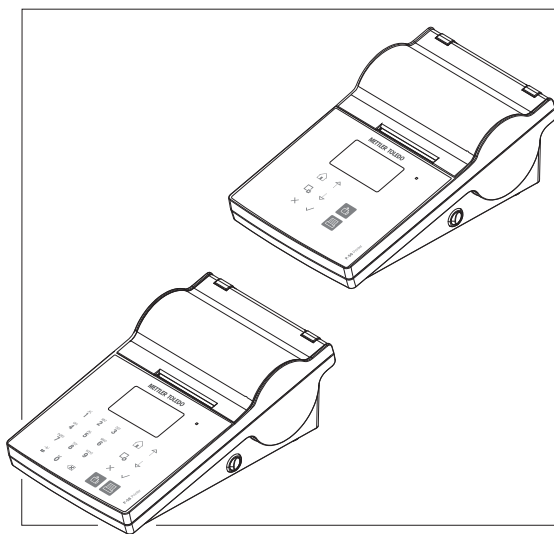
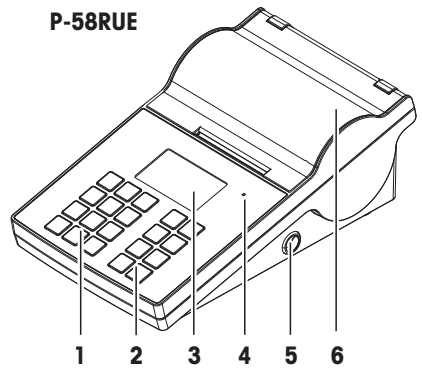
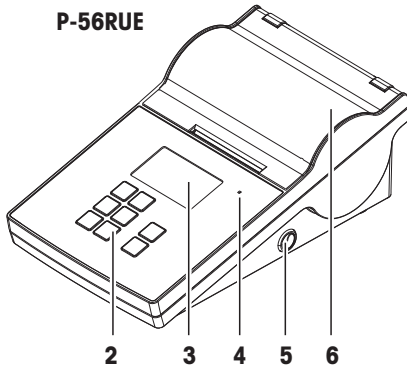
English	User Manual Printer P-56RUE, P-58RUE
Deutsch	Benutzerhandbuch Drucker P-56RUE, P-58RUE
Español	Manual de usuario Impresora P-56RUE, P-58RUE
Français	Guide de l'utilisateur Imprimante P-56RUE, P-58RUE
Italiano	Manuale per l'utente Stampante P-56RUE, P-58RUE
简体中文	用户手册 打印机 P-56RUE, P-58RUE
日本語	ユーザマニュアル プリンタ P-56RUE, P-58RUE
한국어	사용자 매뉴얼 프린터 P-56RUE, P-58RUE

METTLER TOLEDO

Overview

1	Alphanumerical keys	4	Triple status indicator
2	General control keys	5	Power switch
3	Display	6	Printer cover

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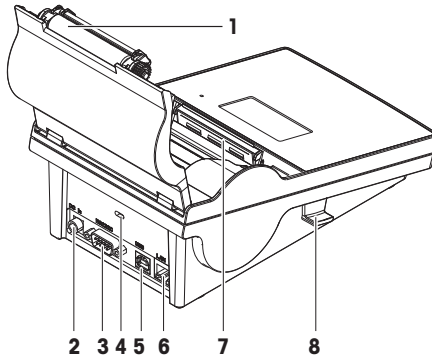
Overview

1	Platen roller	5	USB-B connection
2	Socket for AC/DC adapter	6	Ethernet RJ45 connection
3	RS232 serial interface (9-pin)	7	Thermal printhead
4	Anti-theft slot	8	Handler to open the printer cover

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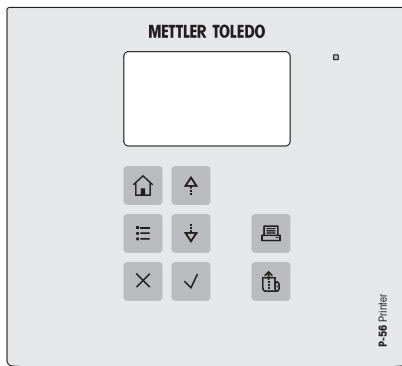
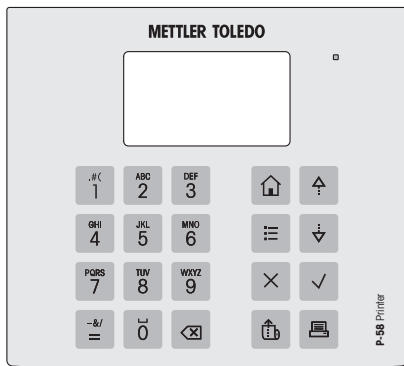
Overview

General Control keys

	Home key	To return to the printer home screen.
	Menu key	To open the section Settings (P-56RUE). To open the section Main menu (P-58RUE).
	Scroll up key	To scroll up through a list or to move the cursor to the left in a string or number entry screen.
	Scroll down key	To scroll down through a list or to move the cursor to the right in a string or number entry screen.
	Cancel key	To cancel a running process/application or the last step.
	Confirmation key	To confirm the current selection.
	Paper feed key	To feed the paper.
	Print key	To print parameters or weighing results.

Alphanumerical keys (P-58RUE)

.#(1	1 . # () , :	PQRS 7	7 P Q R S p q r s
ABC 2	2 A B C a b c	TUV 8	8 T U V t u v
DEF 3	3 D E F d e f	WXYZ 9	9 W X Y Z w x y z
GHI 4	4 G H I g h i	0	0 _
JKL 5	5 J K L j k l	-&/ =	= - & / + * @ % Last entry recall of ID

MNO 6	6 M N O m n o		Delete key To delete the last character.
-----------------	---------------	--	---

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Table of Contents

1	Introduction	3
1.1	Further documents and information	3
1.2	Software download.....	3
2	Safety Information	3
2.1	Definitions of signal words and warning symbols	3
2.2	Product-specific safety information	4
3	Installation and Putting into Operation	5
3.1	Scope of delivery.....	5
3.2	Connecting the printer to the power supply	5
3.3	Inserting the paper roll	5
3.4	Connecting the printer to a METTLER TOLEDO instrument	6
3.4.1	Connecting the printer via the RS232 connection	6
3.4.2	Connecting the printer via the USB connection	7
3.4.3	Connecting the printer via the Ethernet connection.....	7
3.5	Defining language, date and time	7
4	Maintenance	7
5	Technical Data	8
6	Disposal	9

1 Introduction

Thank you for choosing a METTLER TOLEDO accessory. The accessory combines high performance with ease of use.

The P-56RUE and P-58RUE are easy-to-handle direct thermal printers for METTLER TOLEDO instruments. These high-value printers guarantee long-term traceability.

Intended use

The thermal printers P-56RUE and P-58RUE are intended to print data to thermal paper and labels.

1.1 Further documents and information

Finding more information

Search for documents:

▶ www.mt.com/library

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

▶ www.mt.com/contact

1.2 Software download

If needed, download the following software from the METTLER TOLEDO website:

- Ethernet configurator
- Label creator

▶ www.mt.com/printer-software

2 Safety Information

Two documents named "User Manual" and "Reference Manual" are available for this accessory.

- The User Manual is printed and delivered with the accessory.
- The electronic Reference Manual contains the information you need to set up the accessory, put it into operation and maintain it.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.
- The Reference Manual of the main instrument contains additional information on the interface between the accessory and the main instrument.

Only use the accessory according to the User Manual and Reference Manual of the accessory and the Reference Manual of the main instrument. If you do not use the accessory according to these documents or if the accessory is modified, the safety of the main instrument and the accessory may be impaired and Mettler-Toledo GmbH assumes no liability.

2.1 Definitions of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

Signal words

DANGER A hazardous situation with high risk, resulting in death or severe injury if not avoided.

WARNING A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.

CAUTION A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.

NOTICE A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.

Warning symbols

General hazard

Hot surface

Notice

2.2 Product-specific safety information

Your instrument corresponds to the state of the art technology and complies with all recognized safety rules, however, certain hazards may arise in extraneous circumstances. Do not open the housing of the instrument: it does not contain any parts which can be maintained, repaired or replaced by the user. If you ever have problems with your instrument, contact your authorized METTLER TOLEDO dealer or service representative.

Safety notes

WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

WARNING

Injury or damage due to incorrect handling of the battery (placed inside the printer).

- 1 Do not expose the battery or the printer to environmental conditions that exceed those described in the technical data section herein.
- 2 Replace the battery by the same type.
- 3 Dispose the battery or the printer according to the local rules, never damage it or dispose it into fire.

WARNING

Death or serious injury due to electric shock

- 1 Disconnect the printer from the power supply prior to perform maintenance.
- 2 Prevent liquid from entering the printer or AC/DC adapter.

NOTICE

Data Loss

Continuous printing can lead to an overheating of the print head. The printing process stops until the print head has cooled down. During this cooling down phase the printer must not be switched off, otherwise the buffer data will be lost.

NOTICE

Damage to the instrument or malfunction due to the use of unsuitable parts

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

3 Installation and Putting into Operation

Operating conditions

- Use the printer only in dry interior rooms.
For the environmental conditions see the technical data.
- Let the printer acclimate to room temperature.
- Do not use the printer in the presence of humidity and liquids.

3.1 Scope of delivery

The following items are included in the scope of delivery:

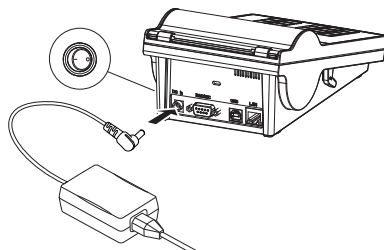
- P-56RUE or P-58RUE thermal printer
- AC/DC adapter
- Country-specific power cable
- USB cable
- RS232 cable
- Set of paper rolls
- User Manual
- Declaration of Conformity

Check the package, the packaging elements and the delivered components for damages. If any components are damaged, please contact your METTLER TOLEDO service representative.

A list of spare parts and accessories can be found in the Reference Manual.

3.2 Connecting the printer to the power supply

- 1 Connect the AC/DC adapter to the socket on the back of the printer and lock it with the screw nut.
- 2 Connect the power cable to the AC/DC adapter and to the power outlet.
- 3 Switch on the printer.

3.3 Inserting the paper roll

CAUTION

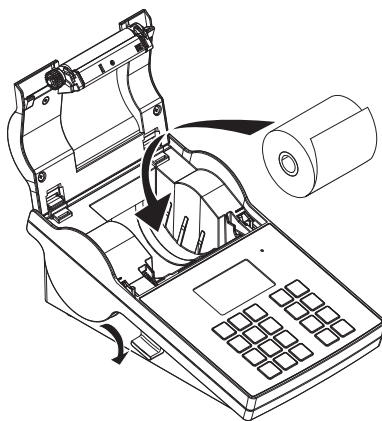
Danger of burns

The thermal printhead is getting hot during the printing process. Wait until the printhead has cooled down before opening the printer cover.


Note

For spare paper rolls, see Accessories and Spare Parts in the Reference Manual.

- 1 Switch on the printer.
- 2 Press down the release key.
➔ The printer cover opens.
- 3 Insert the paper roll into the printer frame.
Note: The paper roll must unroll in the correct direction (see picture).

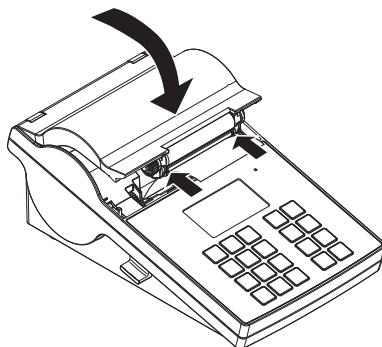- 4 Make sure the paper is in the middle of the paper guides.
- 5 Close the printer cover. Firmly press down until it snaps.
- 6 Press the paper feed key  to check the paper feed.

3.4 Connecting the printer to a METTLER TOLEDO instrument

NOTICE

METTLER TOLEDO recommends connecting the printer to only one instrument at a time. If several instruments are connected METTLER TOLEDO cannot assume the correct functionality of the printer.

3.4.1 Connecting the printer via the RS232 connection

- 1 Connect the printer to the power supply.
- 2 Use the RS232 cable to connect the printer to the METTLER TOLEDO instrument.
Note
The RS232 connection settings of the printer and the instrument must be identical.
- 3 Switch on the printer.

Printer default settings

Baudrate:	9600	Stop bits:	1 Bit
Bits / Parity:	8 / None	Handshake:	XOn/XOff

3.4.2 Connecting the printer via the USB connection

- 1 Connect the printer to the power supply.
- 2 Use the USB cable to connect the printer to a METTLER TOLEDO instrument.
- 3 Switch on the printer.

3.4.3 Connecting the printer via the Ethernet connection

- 1 Connect the printer to the power supply.
- 2 Use an Ethernet cable to connect the printer directly to a METTLER TOLEDO instrument or indirectly over the company's Ethernet.

Note

For installation and configuration of the Ethernet interface you need to have basic knowledge of TCP/IP-based networks, and of network technology in general. If necessary, contact the network administrator of your IT department or your IT support.

3.5 Defining language, date and time

When using the printer for the first time, we recommend to define the printer language, the current time and date.

- 1 Switch on the printer.
 - ➔ The message **Set language & date / time first!** is indicated.
- 2 Use the scroll keys (↓, ↑) to chose a language from the list.
 - ➔ Press the confirmation key ✓.
- 3 Chose a date format.
- 4 Define the date.
- 5 Chose a time format.
- 6 Define the time.
 - ➔ The language, time and date have been defined.

Note

- These settings are retained even if you disconnect your printer from the power supply.
- Date and time will be changed from a connected instrument if the connected instrument supports date and time.
- If the connected instrument supports date and time, the menu items are not available.

4 Maintenance

Cleaning

Make sure to clean the printer regularly.

NOTICE

Damage to the printer

- Do not open the housing of the printer or the AC/DC adapter. They contain no components, which can be cleaned, repaired or replaced by the user.

The printer housing is made of top grade, resistant materials. Therefore, all commercially available mild cleaning agents can be used for cleaning.

Finding more information

See the Reference Manual for further maintenance instructions.

5 Technical Data

Power Supply

AC/DC Adapter:	Primary: 100 V - 240 V, $\pm 10\%$, 50/60 Hz Secondary: 24 V DC, 2.5 A (with electronic overload protection)
Power supply to the printer:	24 V DC, 2.5 A

Protection and Standards

Overvoltage category	II
Degree of pollution	2
Standards for safety and EMC	Declaration of Conformity (part of standard equipment)
Range of application	Indoor use only

Environmental Conditions

Height above mean sea level	up to 4000 m
Ambient temperature range	Operation: 5°C ~ 40°C Storage: -25°C ~ 60°C
Relative air humidity	20% ~ 80% up to 31°C decreasing linearly to 50% at 40°C, non-condensing

Printer Features

Print technology	Direct Thermal
Resolution	8 dots/mm (203 dpi)
Print speed	Max. 101.6 mm/s (4 IPS)
Printable width	54 mm (2.12")
Real time clock	Date and time
Resident fonts	7 Windows bit map fonts: 6 pt, 7 pt, 10 pt, 12 pt, 24 pt, 16x26, OCR A & B
Barcodes	<ul style="list-style-type: none">1-D Bar codes: Code 39, Code 1382-D Bar codes: PDF417 (layout as 1-D), Datamatrix code, QR code, Aztec code
Graphics	<ul style="list-style-type: none">Support BMP and PCXSupport ICO, WMF, JPG, EMF file through softwareSupport image resize, rotating, mapping and inverse through DirectLabel software
Interfaces	<ul style="list-style-type: none">RS232 port (Male DB 9)USB 2.0 device port (B-type)Ethernet 10/100base-TX port (Rj-45)
Control Panel	<ul style="list-style-type: none">Triple color LED: green, orange or red for status indicating8 tactile keys (P-56RUE), 20 tactile keys (P-58RUE)
Display	LCD, backlight, 128 x 64 dots
Display languages	English, German, French, Italian, Spanish, Russian, Chinese (simplified), Japanese, Korean

Materials

Top housing	Die-cast zinc, chromed and plastics (PBT)
Bottom housing	Plastics (PBT)
Platen-roller	

Media

Width	Max. 58 mm (2.28")
Thickness	Min. 0.06 mm (0.003") ~ Max. 0.2 mm (0.008")
Roll diameter	Max. 60 mm (2.36")
Core diameter	12.7 mm (0.5")
Paper roll wound direction	Outwards

USB driver

Link [▶ http://www.fdicchip.com/Drivers/VCP.htm](http://www.fdicchip.com/Drivers/VCP.htm)

Dimension / Weight

Length	220 mm
Height	88 mm
Width	130 mm
Net Weight	1.4 kg, without AC/DC adapter

6 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.

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© 06/2023 METTLER TOLEDO. All rights reserved.