English

## Installation and Operating instructions

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Opzione BTS, 11132535

METTLER TOLEDO

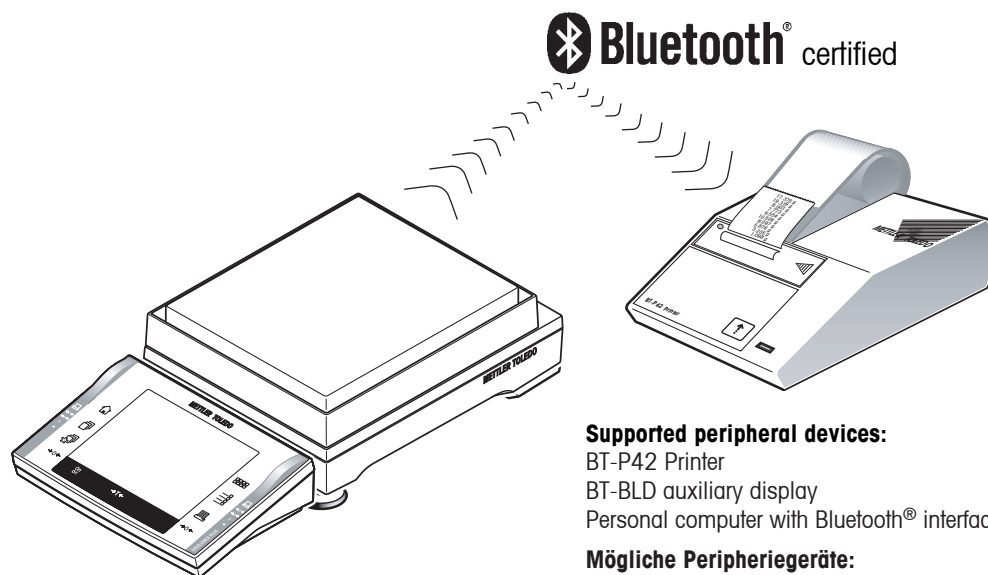
Option BTS is a wireless Bluetooth® interface for connecting a peripheral device to an XP/XS precision balance.

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### Supported peripheral devices:

BT-P42 Printer  
BT-BLD auxiliary display  
Personal computer with Bluetooth® interface

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Numero d'ordine 11132540

Numero d'ordine 11132555

**Technical data:** Wireless data transmission using Bluetooth® technology — Maximum range of 10 m

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Bluetooth is a registered trademark of Bluetooth SIG, Inc.

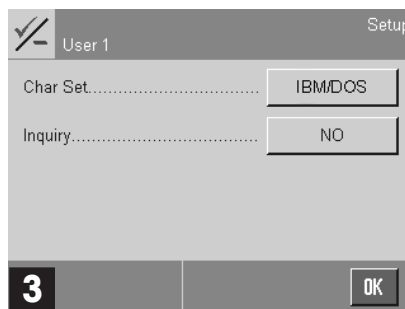
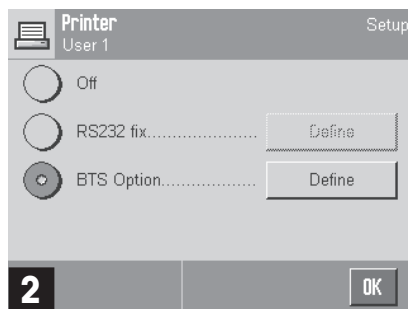
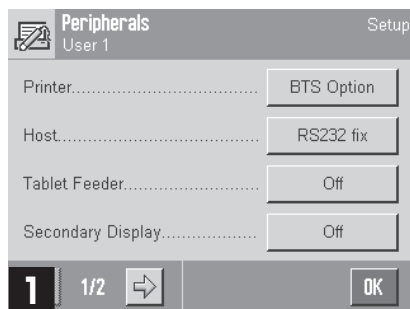
© Mettler-Toledo GmbH 2005 11780627A Printed in Switzerland 0511/6.45

Mettler-Toledo GmbH, Laboratory & Weighing Technologies, CH-8606 Greifensee, Switzerland  
Phone +41-44-944 22 11, Fax +41-44-944 30 60, Internet: <http://www.mt.com>

P11780627

<b>Balance Setup</b>	Menu: System → Peripherie
<b>Waagen-Einstellungen</b>	Menü: System → Peripherie
<b>Réglage de la balance</b>	Menu : Système → Périphériques
<b>Ajustes de la balanza</b>	Menú: Sistema → Periféricos
<b>Impostazioni bilancia</b>	Menu: Sistema → Periferiche

**3 Connect to the peripheral device:** Toggle the [Inquiry] button from "NO" to "YES". Press [OK] to search for and connect to the nearest device. This can take up to 20 seconds. The blue lights on the BTS option of the balance and the peripheral device are continuously lit if the connection is present.

**Important:** Ensure that the only peripheral device within range (maximum of 10 m) of the balance is the one that should be connected to the balance. Devices that are already connected are "busy" and are not detected. They can therefore be positioned within range.

**Note:**

- An existing connection is saved and protected against power failure.
- The connection is interrupted if the interface is changed or the peripheral device is set to "Off". **2**

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