


OptiOx™ Replacement cap
OptiOx™ Messmodul
Module de mesure OptiOx™
Modulo di misurazione OptiOx™
Módulo de medición OptiOx™
OptiOx™ 測量模块
OptiOx™ 測定モジュール
Измерительный модуль OptiOx™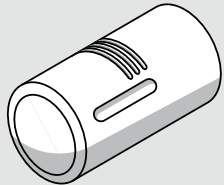

OptiOx™ BOD adapter
OptiOx™ BSB Adapter
Adaptateur DBO OptiOx™
Adattatore BOD OptiOx™
Adaptador BOD OptiOx™
OptiOx™ BOD 适配器
OptiOx™ BOD 測定用アダプタ
Адаптер для определения БПК OptiOx™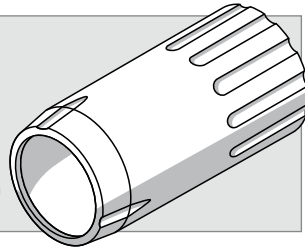

OptiOx™ protective guard
OptiOx™ Schutzaufsatz
Capuchon de protection OptiOx™
Elemento di protezione OptiOx™
Suplemento de protección OptiOx™
OptiOx™ 保护套
OptiOx™ プロテクター
Защитная насадка OptiOx™

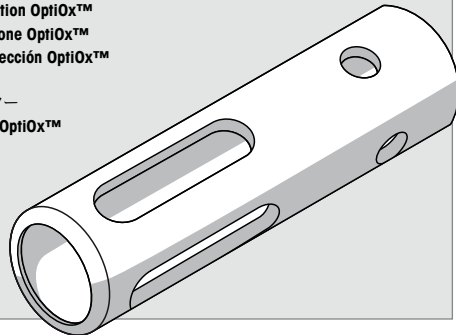GB OptiOx™ sensor equipment

Note: The OptiOx™ sensor contains an internal clock which counts down the 365 day life span of a new sensorcap. The countdown begins once the OptiOx™ cap has been attached, the sensor connected to the device and the first measurement carried out. This process cannot be undone once the first measurement has been carried out.

1. Remove the transport protection from the end of the OptiOx™ sensor. **See A.**
2. Ensure that the two O-rings on the OptiOx™ sensor are correctly positioned. **See B.**
3. Line up the arrow on the OptiOx™ cap with the arrow on the OptiOx™ sensor. **See C.**
4. Push the OptiOx™ cap onto the OptiOx™ sensor until the cap is firmly connected to the sensor. **See D.**
5. Connect OptiOx™ sensor on the mini LTW connection of the device. **See E.**

OptiOx™ sensor storage

Store the sensor in the calibration tube ensuring first that the sponge has been dampened with distilled water or keep the sensor in a beaker filled with distilled water. **See F.**

Note: Detailed information on the OptiOx™ sensor can be found in the user manual for the device.

 This product must not be disposed of with domestic refuse. Please dispose of this product in accordance with local regulations at a separate collection point for electrical and electronic equipment.

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