
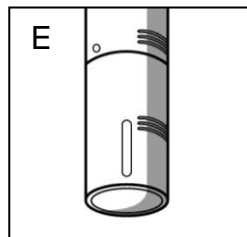
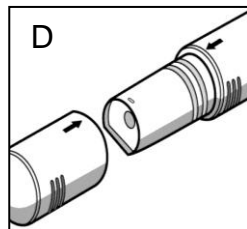
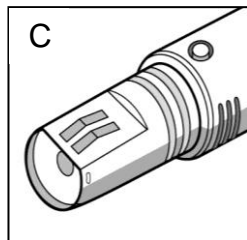
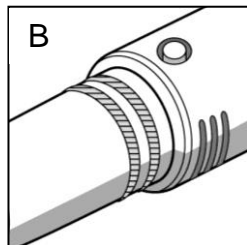
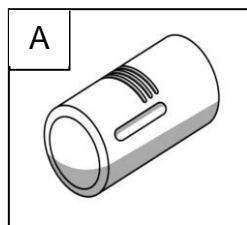
**OptiOx™ Replacement Cap
(P/N 51343164)**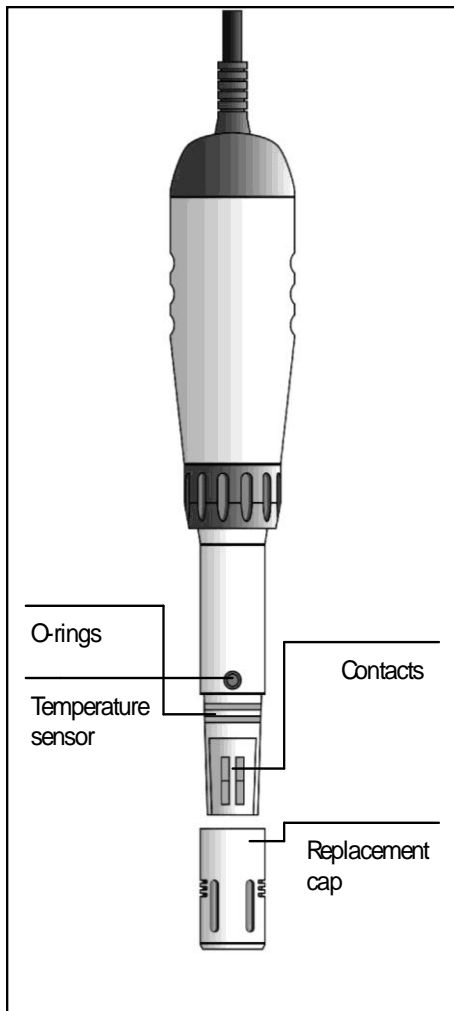

Replacement of the OptiOx™ sensor cap

The InLab® OptiOx sensor contains an internal clock which counts down the 365 day life span of a new sensor cap. The countdown begins once the OptiOx™ cap has been attached, the sensor connected to the device and the first measurement carried out. This process cannot be undone once the first measurement has been carried out.

Note: Install the cap by the date printed on the package. Keep the cap in its sealed packaging until you are ready to install it. Install promptly.

1. Pull the cap off the sensor without twisting (A).
2. Remove the existing O-rings from the sensor (B).
3. Use a lint-free cloth to remove any moisture from the sensor body. Make sure that there is no moisture in the O-ring grooves (C).
4. Use your finger to apply a layer of lubricant around the O-ring grooves. Place the O-rings on the sensor. Apply another thin layer of lubricant to the O-rings and grooves.

Do not transfer grease to the lens or contacts!

5. Gently wipe the lens with a clean lens cloth and allow it to dry thoroughly. Examine the lens for scratches or dirt.
6. Remove the new cap from its sealed packaging. Line up the arrow on the cap with the arrow on the sensor (D).
7. Push the cap onto the sensor until it is firmly connected. Make sure that the O-rings are not pinched or rolled between the cap and sensor (E).
8. After replacing the cap, a water-saturated air calibration should be performed.

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