

MA Balance

Software Release Notes

Contents

- 1 Software Version History 1
- 2 Updating the Balance Software 2
- 3 Software Release Information 3
- 4 Support..... 3

1 Software Version History

Date	Version	Topic	Description	New TDNR
Dec. 11, 2023	1.1.0	Initial release	Read more in SW information V 1.1.0	-

2 Updating the Balance Software


Search for software:

▶ <http://www.mt.com/labweighing-software-download>

Contact a local METTLER TOLEDO service representative if you need support updating the software.

Software update process

The software update is only accessible to users with the corresponding rights. Make sure that the USB storage device used contains only one software in the file format MOT.

- A USB storage device containing the software is connected to the balance and not to be removed during the software update.
- 1. Press  to enter the settings menu.
- 2. Navigate to the menu entry **BAL.SET**.
- 3. Navigate to the menu entry **MAINT**.
- 4. Navigate to the menu entry **SOFTWARE UPDATE**.
 - ➔ A request to insert a USB storage device is displayed.
- 5. Insert a USB storage device with the desired software version.
- 6. Select **START** to run the update.
- ➔ The balance performs the update. After completing the update, the balance reboots.

3 Software Release Information

SW Version 1.1.0

Date	Version	Update History	New TDNR
Dec. 11, 2023	1.1.0	New	Initial version for MA

4 Support

For any further support, please contact us: www.mt.com/contact