InFit 76X

Stationary Housings

Quick Setup Guide

Contents

cs	Průvodce rychlým nastavením	2
da	Lynvejledning	5
de	Quick-Setup-Leitfaden	8
en	Quick Setup Guide	11
es	Guía de configuración rápida	14
fi	Pika-asetusopas	17
fr	Guide de paramétrage rapide	20
hu	Gyorsbeállítási útmutató	23
it	Guida alla configurazione rapida	26
ja	クイックセットアップガイド	29
ko	빠른 설정 가이드	32
nl	Beknopte handleiding	35
pl	Podręcznik szybkiej konfiguracji	38
pt	Guia de configuração rápida	41
ru	Руководство по быстрой настройке	44
sv	Snabbinstallationsguide	47
th	คู่มือการตั้งค่าอย่างรวดเร็ว	50
zh	快速设置指南	53

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Quick Setup Guide for InFit 76X

1 Introduction

The METTER TOLEDO InFit™ 76X stationary housings are intended for use in conjunction with specified METTLER TOLEDO electrodes/sensors, namely pH and redox (ORP) combination electrodes as well as dissolved oxygen (DO), CO₂, conductivity or turbidity sensors. Please read through this quick set-up guide and the operating instructions carefully before commissioning, in order to ensure trouble-free use. Operating and maintenance should be carried out only by trained personnel and staff who have read and understood the operating instructions.

2 Safety Instructions

 **Notice:** Before every start-up, the housing must be checked for:

- Damage to the connections, fastenings, etc.;
- Leakage;
- Perfect functioning;
- Authorization for use in conjunction with other plant equipment and resources.

 **Notice:** The manufacturer / supplier accepts no responsibility for any damage caused by unauthorized attachments or for the incorporation of spare parts which are not of METTLER TOLEDO provenance. The risk is borne entirely by the operator.

 **Caution:** Before the housing is put into operation, the operator must have already clarified that use of the sensor in conjunction with the other associated equipment and resources is fully authorized.

 **Caution:** A defective housing may neither be installed nor put into operation. Faulty containment or installation out of conformance with regulations and instructions can lead to the escape of medium or to pressure surges (explosion), potentially harmful both to persons and to the environment.

 **Caution:** If an electrode/sensor is broken it must be replaced immediately.

 **Warning:** Before dismantling the housing or commencing any maintenance work on it, ensure that the equipment in which the housing is installed is in a safe condition (depressurize, no explosion risk, empty, rinse, vent, etc.).

 **Warning:** Please observe, read and fully understand the special conditions of use for installation in Ex classified areas.

3 Product Description

The label on the housing contains the following information:

METTLER TOLEDO www.mt.com Im Hackacker 15, CH-8902 Urdorf, Switzerland	Manufacturer of the housing and contact details
P/N	Part number
S/N	Serial number
PS	Maximum permissible pressure
TS	Maximum permissible temperature
⊕ II 1/2G Ex h IIC T6...T3 Ga/Gb ⊕ II 1/2D Ex h IIIC T69°C...T141°C Ga/Gb	ATEX tested and description (model dependent)
CE _{xxxx}	Ex classification testing notified body number
See instruction manual	See instruction manual for detailed description

4 Installation and Commissioning

1. While unpacking, check the housing for mechanical damage. Please immediately report any damage to your local METTLER TOLEDO representative.
2. Isolate the pipe/tank/vessel etc. from the housing (depressurize, empty, rinse, purge and vent, etc.) before starting any installation work.
3. Fit the housing to flange/NPT thread of process connection.
4. Fit the electrode/sensor to the housing and connect the electrode/sensor and the transmitter with corresponding cable.

 **Note:** Perform step #4 followed by #3 for InFit 762/763 housing.

5. Adjust compensation pressure: The compensation pressure may be adjusted at the valve insert of the pressure gauge using the supplied air pump, or by connecting a permanent, oil-free and filtered pressure supply (for housing InFit 763 or InFit 764 only).

Refer to respective housing instruction manual for detail description.

5 Operation: Inspection Work

1. Visually check fastenings (ring nut, flange, NPT thread) of the housing at the process connection for firm sealing and leaks.
2. If necessary, check the condition of the electrode/sensor (refer to sensor instruction manual for cleaning and calibration procedures).
3. Calibrate/clean the electrode/sensor if necessary (refer to sensor and transmitter instruction manuals).

6 Maintenance

1. Isolate the pipe/tank/vessel etc. from the housing (depressurize, empty, rinse, purge and vent, etc.).
2. Remove the sensor from the housing.
3. Remove the stationary housing from the pipe/tank/vessel etc. (if changing of housing seal is needed).

 **Note:** Perform step #3 followed by #2 for InFit 762/763.

4. Change the housing seal.
5. Calibrate/clean the electrode/sensor with reference to sensor and transmitter instruction manuals.
6. Follow the installation and commissioning procedure.

7 Environmental protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

8 Warranty

On manufacturing defects, 12 months after delivery.

9 Storage Condition

We recommended following the shutdown procedure as described in the housing instruction manual. The cleaned and dried housing should be stored at room temperature.

10 EC Declaration of Conformity

The complete declaration is available in the declaration of conformity certificate.

InFit is a trademark of the METTLER TOLEDO Group.

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All rights reserved