# Quick Setup Guide Transmitter M100/2H DR

---

## Multilingual Quick Setup Guide

cs	Průvodce rychlým nastavením	3
da	Lynvejledning	13
de	Quick-Setup-Leitfaden	23
en	Quick Setup Guide	33
es	Guía de configuración rápida	43
fi	Pika-asetusopas	53
fr	Guide de paramétrage rapide	63
hu	Gyorsbeállítási útmutató	73
it	Guida alla configurazione rapida	83
ja	クイック セットアップ ガイド	93
ko	빠른 설정 가이드	103
nl	Beknopte handleiding	113
pl	Podręcznik szybkiej konfiguracji	123
pt	Guia de Configuração Rápida	133
ru	Руководство по быстрой настройке	143
sv	Snabbinstallationsguide	153
th	คู่มือการตั้งค่าอย่างรวดเร็ว	163
zh	快速操作指南	173

**METTLER TOLEDO**

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---

### Obsah

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---

### Content

1	Safety and Intended Use	34
2	Technical Data and Further Information	35
3	Mounting	35
4	Electrical Connection	35
5	Menu Structure	37
6	General Setup	38
7	Sensor Calibration	41
8	Maintenance	42
9	Disposal	42
10	EC-Declaration of Conformity	42

# 1 Safety and Intended Use

 **Note:** The Quick Setup Guide is a brief operating instruction.

The M100 transmitter must only be installed, connected, commissioned, and maintained by qualified specialists e.g. electrical technicians in full compliance with the instructions in this Quick Setup Guide, the applicable norms and legal regulations.

The specialist must have read and understood this Quick Setup Guide and must follow the instructions it contains. If you are unclear on anything in this Quick Setup Guide, you must read the Operating Instructions (supplied on CD-ROM). The Operating Instructions provides detailed information on the instrument.

The M100 transmitter should be operated only by personnel familiar with the transmitter and who are qualified for such work.

## Intended Use

The M100 DIN rail transmitter is a 2-wire transmitter for analytical measurements with HART communication capabilities. The M100 is a 1-channel, multi-parameter unit for pH/ORP, conductivity and dissolved oxygen measurement. It is only compatible with ISM® sensors.

The M100 transmitter is designed for use in the process industries, in non-hazardous areas.

## M100 parameter fit guide

Parameter	M100/2H DR
	ISM
pH/ORP	•
pH/pNa	•
Conductivity 4-e	•
Amp. Dissolved Oxygen ppm / ppb / trace	• <sup>1)</sup> / • <sup>2)</sup> / • <sup>2)</sup>

1) INGOLD and THORNTON sensors

2) THORNTON sensors

## 2 Technical Data and Further Information

The most important technical data such as supply voltage are given on the name plate on the transmitter. For further technical data such as accuracy refer to the Operating Instructions. This document, the Operating Instructions and software are on the supplied CD-ROM. You can also download the documentation via Internet "[www.mt.com/M100](http://www.mt.com/M100)".

## 3 Mounting

The M100 DIN rail transmitter is suitable for 35 mm wide DIN rail systems. For installation drawings refer to the Operating Instructions.

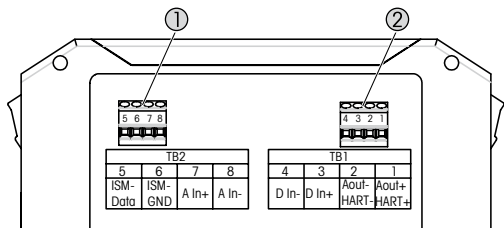
## 4 Electrical Connection

 **Note:** Power off instrument during electrical connection.

1. Switch off supply voltage.
2. Connect mains supply (14 to 30 V DC) to the terminals **1** and **2**.
3. Connect sensor to terminals **5** and **6**.
4. Connect analog output and HART signal to terminals **1** and **2**.
5. Connect digital input for switching transmitter in Hold state to terminals **3** and **4**.
6. Connect analog input signal 4 to 20 mA for pressure compensation to terminals **7** and **8**.


For terminal definitions see next page.

## Terminal Definition

- 1 Terminal block TB2
- 2 Terminal block TB1

## LED Functionality


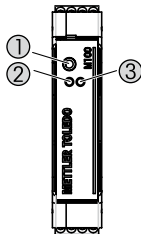
- 1 Port: Interface for service  
e.g. firmware update
- 2 LED green
- 3 LED red

Terminal			Definition	
TB2	5	ISM-Data	Sensor input, see "Safety and Intended Use" chapter on Page 34.	
	6	ISM-GND		
	7	Ain+		Analog input 4 to 20 mA (for pressure compensation)
	8	Ain-		
TB1	4	DI-	Digital input (for switching transmitter in Hold state)	
	3	DI+		
	2	Aout-, HART-	Power connection 14 to 30 V DC, Analog output and HART signal	
	1	Aout+, HART+		

LED Green	LED Red	Definition
X	–	Operation correct
X	X	Error
X	Blinking: 0.5 s: on 0.5 s: off	Warning

## 5 Menu Structure
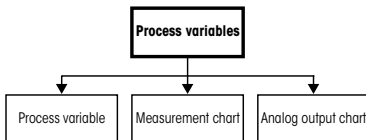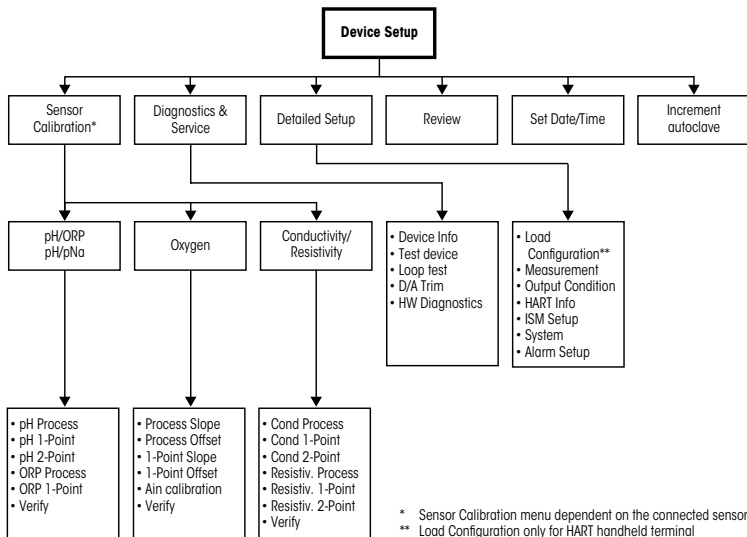
## 6 General Setup

You configure the M100 transmitter either via a configuration tool, an asset-management-tool or via a HART handheld terminal.

Prerequisite: The M100 transmitter and the sensor are mounted and electrically connected.

### Configuration tool or asset-management tool

**Note:** The DTM and the configuration tool PACTWare™ are on the supplied CD-ROM. You can also download the DTM via Internet “[www.mt.com/M100](http://www.mt.com/M100)”.

For steps 1 to 5 and step 13 refer to the documentation of the tool.

1. Install configuration tool e.g. PACTWare™ or asset-management-tool.
2. Install DTM for HART interface and DTM for M100 transmitter.
3. Update device catalog.
4. Build up connection. Check COM port settings if necessary.
5. Select **Sensor Type**.  
Menu path: Detailed Setup > Measurements > Channel Setup > Sensor Setup.
6. Load configuration from device.
7. Set **Tag** or/and **Long Tag**.  
Menu path: Device Setup > Detailed Setup > HART Info
8. Set **Date** and **Time**. Set the time in the 24 hour format. The time format cannot be changed. Menu path: Device Setup > Set Date/Time
9. Set range of the analog output signal.  
Menu path: Detailed setup > Output Condition > Analog output > Range
  - **URV** (Upper Range Value) and **LRV** (Lower Range Value): The values can be changed to the current measurement range. The values have to be within the measurement limits of the sensor.
  - **USL** (Upper Sensor Limit) and **LSL** (Lower Sensor Limit): The limits are defined by the sensor and cannot be changed.

10. Define the process variables **PV**, **SV**, **TV** and **QV**  
Menu path: Device Setup > Detailed Setup > Measurements > Channel Setup
11. Calibrate sensor. Menu path: Device Setup > Sensor Calibration
12. Perform further settings. See Operating Instructions of the M100 transmitter.
13. Store configuration to device.

## HART handheld terminal

**Note:** The DD "008E8E7F0101.hhd" is on the supplied CD-ROM. You can also download the DD via Internet "[www.mt.com/M100](http://www.mt.com/M100)".

For "Step 1" refer to the documentation of the HART handheld terminal.

1. Check if the DD of the M100 transmitter has already been installed on the HART handheld terminal. Install the DD if necessary.
2. The communication is built up automatically.
3. Select **Sensor Type**.  
Menu path: Detailed Setup > Measurements > Channel Setup > Sensor Setup.
4. Load configuration from device. Menu path: Device Setup > Detailed Setup
5. Set **Tag** or/and **Long Tag**.  
Menu path: Device Setup > Detailed Setup > HART Info
6. Set **Date** and **Time**. Set the time in the 24 hour format. The time format cannot be changed. Menu path: Device Setup > Set Date/Time
7. Set range of the analog output signal.  
Menu path: Detailed setup > Output Condition > Analog output > Range
  - **URV** (Upper Range Value) and **LRV** (Lower Range Value): The values can be changed to the current measurement range. The values have to be within the measurement limits of the sensor.
  - **USL** (Upper Sensor Limit) and **LSL** (Lower Sensor Limit): The limits are defined by the sensor and cannot be changed.
8. Define the process variables **PV**, **SV**, **TV** and **QV**  
Menu path: Device Setup > Detailed Setup > Measurements > Channel Setup

9. Calibrate sensor. Menu path: Device Setup > Sensor Calibration
10. Perform further settings. See Operating Instructions of the M100 transmitter.

### **Change Sensor Type**

Prerequisite: The M100 transmitter and another sensor type are mounted and electrically connected.

If you change the sensor type e.g. a Conductivity sensor with a pH sensor, perform the following procedure:

1. Built up connection.
2. Select **Sensor Setup**.  
Menu path: Detailed Setup > Measurements > Channel Setup > Sensor Setup.
3. Start "Sensor Setup".
4. Select the new sensor type for the **Sensor Type** parameter.
5. Load configuration from device.
6. The configuration in the configuration tool or in the HART handheld is updated. If you have selected the correct sensor type, the "Verify" menu will be shown.

## 7 Sensor Calibration

(PATH: Device setup/Sensor Calibration)

**Note:** For best process calibration results observe the following points. Take grab sample as close as possible to the measurement point of the sensor. Measure the sample at process temperature.

Via configuration tool, asset-management-tool or HART handheld terminal you can calibrate the sensor with the “Process”, “1-Point” or “2-Point” calibration method. For the calibration methods see Operating Instructions of the M100 transmitter. Via iSense software you can calibrate the sensor with the “1-Point” or “2-Point” calibrate method. See Operating Instructions of the iSense software.

As soon as the calibration is in progress no other calibration can be started.

### Sensor Calibration Menu

For the calibration methods see “Menu Structure” chapter on Page 37.

After every successful calibration the following options are available:

- **Adjust**

Calibration values are adopted and used for the measurement. Additionally, the calibration values are stored in the calibration history.

- **Calibrate**

Calibration values are stored in the calibration history for documentation, but not be used for the measurement. The calibration values from the last valid adjustment are further used for the measurement.

- **Abort**

Calibration values are discarded.

## **8 Maintenance**

The transmitter requires no maintenance.

Clean the surfaces with a soft damp cloth and dry the surfaces with a cloth carefully.

## **9 Disposal**

Observe the applicable local or national regulations concerning the disposal of "Waste electrical and electronic equipment".

Disassemble the transmitter according to resources. Sort resources and supply them to recycling. Non-recyclable materials have to be disposed of in an environmental-friendly manner.

## **10 EC-Declaration of Conformity**

The EC-Declaration of conformity is part of the delivery.

ISM is a registered trademark of the Mettler-Toledo Group in Switzerland, Brazil, USA, China, European Union, South Korea, Russia and Singapore.

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© Mettler-Toledo GmbH 08/2016.
- D** Mettler-Toledo GmbH, Prozeßanalytik, Ockerweg 3, D-35396 Gießen
- E** Mettler-Toledo S.A.E., C / Miguel Hernández, 69-71, ES-08908 L'Hospitalet de Llobregat (Barcelona)
- F** Mettler-Toledo, Analyse Industrielle S.A.S., 30, Boulevard Douaumont, F-75017 Paris
- CN** Mettler-Toledo International Trading (Shanghai) Co. Ltd., 589 Gui Ping Road, Cao He Jing
- UK** Mettler-Toledo LTD, 64 Boston Road, Beaumont Leys, GB-Leicester LE4 1AW