# InSUS 607

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---

## 1 Introduction

Thank you for buying this METTLER TOLEDO product. InSUS™ optical dissolved oxygen (DO) sensors are gamma irradiation-sterilizable single-use DO sensors. The sensors have been designed especially for deployment in single-use bags used in biotechnical processes. The sensors offer a design that allows safe and easy installation in a standard Eldon James port disc with a 1" barb. This port disc is welded into the bag film prior to sensor installation.

## 2 Safety Instructions

The sensor should be handled carefully. Local safety regulations are applicable to the handling of the sensor. Installation and operation should be carried out only by trained personnel and staff familiar with the unit and who have read and understood this installation manual.

## 3 Product Description

InSUS 607

- 1** Protective cap on M16x1 thread  
(for protection of inner surface of optical spot from ambient light during sensor transport and storage)
- 2** Retainer ring
- 3** O-rings (EPDM)
- 4** Sensor shaft (HDPE) with DO sensing optical spot assembled on sensor tip

## 4 Installation

### 1. Unpacking

On unpacking, check the sensor for possible mechanical damage of the O-rings and the M16x1 thread. Report any damage to your METTLER TOLEDO supplier.

### 2. Installation into an Eldon James bag port

**NOTICE:** 1" Eldon James ports are available in different polymer materials. For a leak proof sensor installation METTLER TOLEDO offers InSUS sensors with different O-ring configurations. If in doubt of the correct configuration for your requirements, please contact your METTLER TOLEDO supplier.

Recommended port/sensor combinations are:

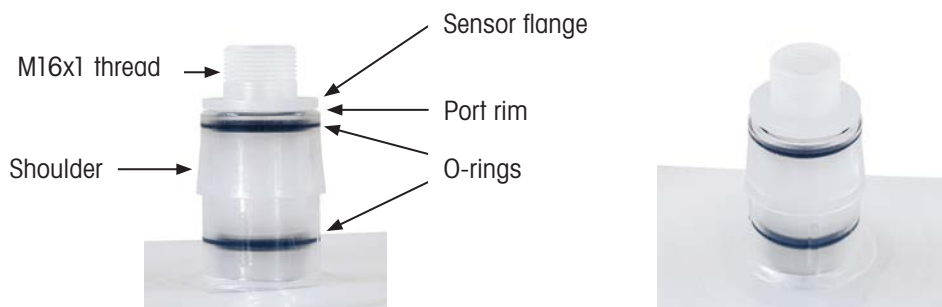
1" Eldon James port material	Compatible InSUS sensor types
Engage™ 8402	InSUS 607
Dowlex™ 2517	InSUS 607/S

If applicable in your production environment use a mixture of 70 % isopropyl alcohol and 30 % deionized water for carefully lubricating the sensor O-rings prior to insertion into the Eldon James port.

Remove the black protective cap and the retainer ring from the sensor. Push the sensor all the way into the port until the sensor flange is flush with the port rim. This is ideally done with a slight twisting motion to facilitate the insertion.


**NOTICE:** Make sure that the O-rings are still located completely in their grooves after this insertion procedure. If this is not the case, pull the sensor out of the port, reposition the O-rings and repeat the insertion procedure.

Correct sensor and O-ring positions after sensor insertion into an Eldon James port.

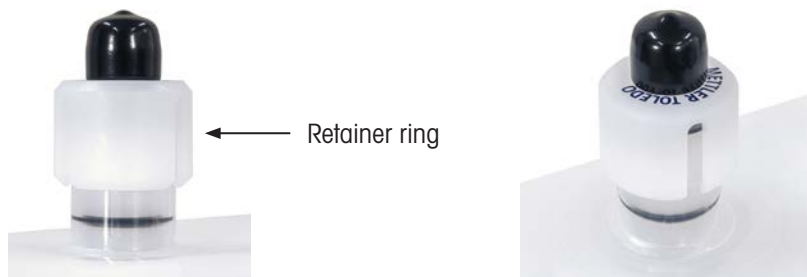Push the retainer ring over the sensor flange and port shoulder until it **clicks** into place.

Correct position of the retainer ring and the black protective cap on sensor and bag port.

**CAUTION:** Ensure that the black protective cap is installed back on the sensor thread for bag storage and shipment!

## 5 Environmental Protection

Please recycle where facilities exist. Check with your Local Authority for disposal or recycling advice.

## 6 Warranty

In the event of manufacturing faults, 12 months warranty from date of production is granted.

InSUS is a trademark of the METTLER TOLEDO Group.

Engage and Dowlex are trademarks of the Dow Chemical Company.

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All rights reserved