# METTLER TOLEDO DG111-SC

**Glass electrode**  
**Glaselektrode**  
**Electrode de verre**  
**Electrodo de vidrio**  
**Стекланный электрод**  
**玻璃电极**

## **Combined pH electrode**

with ceramic frit for direct pH measurements and acid/base titrations in aqueous solutions

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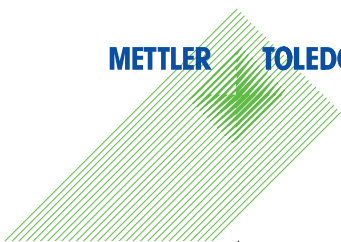
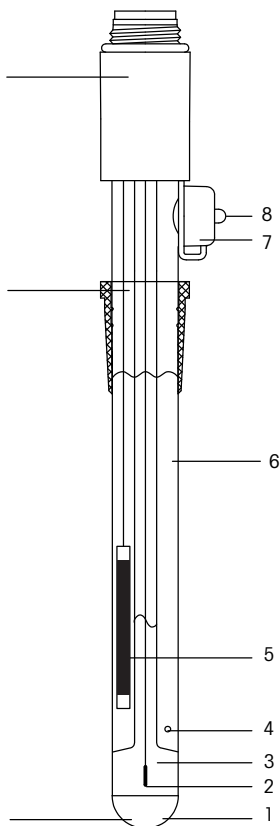
## 1. Pre-operation treatment

- 1.1 The internal buffer solution (3) must cover the whole inside of the glass membrane (1). Remove air bubbles by gently shaking in vertical direction.
- 1.2 To equalize pressure remove the rubber cap (7) or perforate its nipple (8).
- 1.3 Refill the reference electrolyte (6) – (e.g. METTLER TOLEDO No. 51340049) – to approx. 1 cm below the aperture (7).
- 1.4 To activate the glass membrane place the electrode in a tenfold diluted reference electrolyte solution. Use a 0.5 mol/L KCl solution if no reference electrolyte is at hand.

## 2. Adjustment and calibration

The parameters zero-point and slope are measured by adjustment. A new electrode should have parameter values according to 2.1–2.2. Instructions for the adjustment method are given in the titratormanual.

- 2.1 The zero point of the electrode (pH value at  $E = 0$  mV) is approx. pH 7. The electrode potential, measured in a buffer solution of pH 7, must be between  $-30$  mV and  $+30$  mV.
- 2.2 The electrode slope (potential change per pH unit) at the pH range 4–9 must be greater than  $-55$  mV/pH (at 25 °C).
- 2.3 You should test the response time of the electrode. After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 2.4 If calibrating a used electrode the zero point should be within pH 6 and 8, the slope not less than  $-52$  mV/pH, and the potential change after 60 s should not exceed 3 mV/30 s.

## 3. Storage and lifetime

- 3.1 Store the electrode in the reference electrolyte – the frit (4) has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 3.2 Never let the glass membrane dry out, otherwise it has to be reactivated and the KCl crystals in/on the frit have to be dissolved.
- 3.3 The lifetime of the electrode is 6 months to 3 years, depending on its use and treatment.

#### 4. Maintenance instructions

- 4.1 The measuring range extends from pH 0 to 14, the tolerable temperature range from 0 to 80 °C (32–176 °F).
- 4.2 The sample solution should never enter through the ceramic frit into the reference electrode. Therefore the level of the reference electrolyte must always be higher than the sample solution level.
- 4.3 No reaction should take place between the sample solution and the reference electrolyte: sulfides, bromides, iodides and cyanides react with Ag<sup>+</sup> and some cations (Ag<sup>+</sup>, Hg<sup>2+</sup>, Au<sup>3+</sup>, Pb<sup>2+</sup>) react with Cl<sup>-</sup> of the reference electrolyte and cause a blockage of the frit.
- 4.4 Solutions containing protein quickly block the frit, due to protein precipitates.
- 4.5 Just blot the electrode after rinsing, never rub it dry.
- 4.6 Do not let the cable connector plug corrode.
- 4.7 A dried out glass membrane can be reactivated by placing it for several hours in 0.1 mol/L HCl.

#### 5. Cleaning procedures

- 5.1 Contamination of the reference electrolyte: Drain the reference electrolyte through the aperture (7) and proceed according to 1.3 and 1.4.
- 5.2 Frit blocked by AgCl: place it overnight in concentrated ammonia, rinse and condition it for 1 hour in a buffer solution pH 4.
- 5.3 Frit blocked by sulfide precipitate: place it in a solution of thiourea (8% thiourea/1 mol/L HCl, e.g. METTLER TOLEDO No. 51340070) until the dark color has vanished.
- 5.4 Frit blocked by protein: place it at least for 1 hour in pepsin solution (5% pepsin/0.1 mol/L HCl, e.g. METTLER TOLEDO No. 51340068).
- 5.5 Other contaminations: place it for a few minutes in an ultrasonic bath with deion. water, ethanol or a mixed acid solution (e.g. aqua regia).

#### 6. Regeneration

If the electrode does not meet the qualifications according to 2.4 anymore, try a corrosion of the glass membrane: place it 1–2 minutes in diluted HF solution (2% HF/5% HCl, e.g. METTLER TOLEDO No. 51340073), rinse thoroughly with deion. water and treat it as in 5.1.

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