Jewelry JL-G
Jewelry JL-G/L

English

Compact electromechanical balances determine the weight (mass) by measuring the exerted gravity force of the weighing good. As the gravity factor depends on latitude and altitude, the balance must necessarily be adjusted at the location of use. This is done with the delivered external reference weight, according to this installation instruction.

If the balance is used for trade (price based on weighing result) the balance must be switched to the approved mode. A Weights and measures representative inspects and secures the balance against manipulation and fraud, means securing stickers or/and wire seal.

To avoid any conflict with the legislator, every end-user is requested to perform the above steps. Depending on state or country legislation, omission is severely punished.

See procedure described on the next page.

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**ATTENTION**

Service mode can only be selected when the balance is operated with the AC adapter.

**Step 1**

- Press and hold to switch off the balance.
- Press calibration switch **C** until **SERVICE** is displayed (e.g. using a paperclip).

If no key is pressed for 30 minutes, the balance automatically reverts from service mode back to weighing mode.

Step 2

- Press and hold until **METROL** is displayed.
(Press to abort or exit.)

Step 3

- Choose **noAPPr** with and save it by pressing and holding .

Step 4

- Press and hold until **MENU** is displayed.
 - Change the value to the supplied adjustment weight by pressing and save it by pressing and holding .
- (Press to abort or exit.)

Step 5

- Switch off and switch on by pressing .
- Adjust (calibrate) the balance by pressing and holding and follow the instructions on the display.
- Set the balance to the approved mode by repeating steps 1 and 2.
- Repeat Step 3, choose **APPr** and save it by pressing and holding .

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