Quick Setup Guide

Transmitter M400

Multilingual Quick Setup Guide

English	3
Deutsch	19
Français	35
Italiano	51
Español	67
Português	83
Русский	99
日本語	115

METTLER TOLEDO

A graphic element consisting of a series of parallel diagonal lines that form a stylized arrow pointing downwards and to the right, positioned below the company name.

Quick Setup Guide

Transmitter M400

Content

1	Operation	4
2	Menu Structure	5
3	Terminal Block (TB) definitions	6
4	Wiring example for pH Transmitter	8
5	General Setup (applies for all parameters)	9
6	pH Calibration incl. ISFET	10
7	O ₂ Calibration	12
8	CO ₂ Calibration	13
9	M400 Type 1 Cond Ind only: Conductivity Calibration for inductive Sensors	15
10	Conductivity Quick Setup	17
11	Conductivity Calibration for 2-e and 4-e Sensors	17

1 Operation

Entry of data values, selection of data entry options

Use the ▲ key to increase or the ▼ key to decrease a digit. Use the same keys to navigate within a selection of values or options of a data entry field.

Note: Some screens require configuring multiple values via the same data field (ex: configuring multiple setpoints). Be sure to use the ► or ◀ key to return to the primary field and the ▲ or ▼ key to toggle between all configuration options before entering to the next display screen.

Navigation with ↑ in Display

If a ↑ is displayed on the bottom right hand corner of the display, you can use the ► or ◀ key to navigate to it. If you click [Enter] you will navigate backwards through the menu (go back one screen). This can be a very useful option to move back up the menu tree without having to exit into the measuring mode and re-enter the menu.

Exit menu

Note: Exit the menu at any time by pressing the ◀ and ► key simultaneously (ESCAPE). The transmitter returns to the Measurement mode.

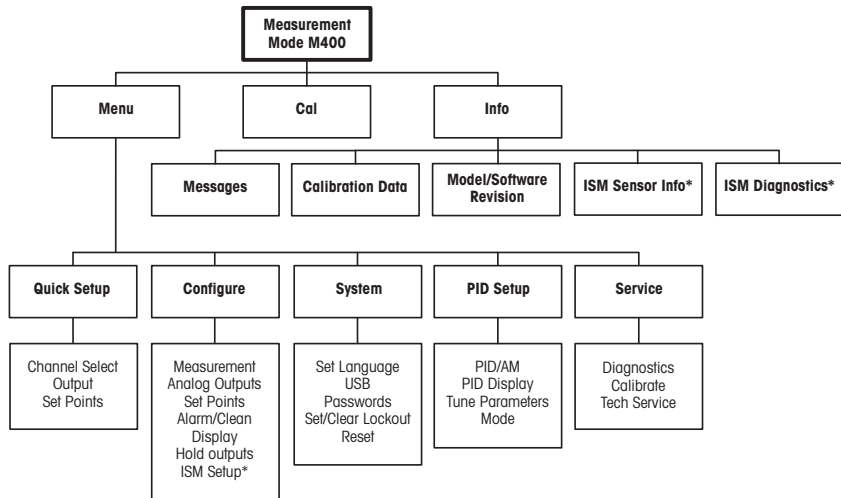
“Save changes” dialog

Three options are possible for the “Save changes” dialog:

- “Yes & Exit”: Save changes and exit to measuring mode
- “Yes & ↑”: Save changes and go back one screen
- “No & Exit”: Don’t save changes and exit to measuring mode

The “Yes & ↑” option is very useful if you want to continue configuring without having to re-enter the menu.

2 Menu Structure


* Only available in combination with ISM sensors.

3 Terminal Block (TB) definitions

Power connections are labeled –N for Neutral and +L for Line, for 100 to 240 VAC or 20–30 VDC.

TB2 for ½ DIN	
1	AO1+
2	AO1–/AO2–
3	AO2+
4	AO3+
5	AO3–/AO4–
6	AO4+
7	DI1+
8	DI1–/DI2–
9	DI2+

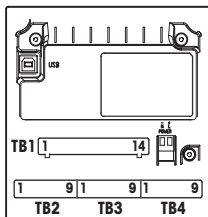Note: This is a 4-wire-product with an active 4–20 mA analog output. Please do not supply to Pin1–Pin6 of TB2.

TB1 for ½ DIN			
1	NO1	8	NC5
2	COM1	9	COM6
3	NC1	10	NO6
4	NO2	11	NO3
5	COM2	12	COM3
6	NC2	13	NO4
7	COM5	14	COM4

NO = normally open (contact is open if unactuated).

NC = normally closed (contact is closed if unactuated).

TB4 – ISM (digital) Sensors

Term.	Function	Optical Oxygen Color	pH, amp. Oxygen, Cond 4-e Color
1	24 VDC	brown	–
2	GND (24 VDC)	black	–
3	1-Wire	–	transparent (cable core)
4	GND (5 VDC)	grey and yellow	red (shield)
5	–	–	–
6	GND (5 VDC)	–	–
7	RS485-B	blue	–
8	RS485-A	white	–
9	5 VDC	–	–

TB3 – Conventional (analog) Sensors

Term.	Cond 2-E or 4-E		O ₂	
	Function	Color*	Function	Color
1	Cnd inner1	white	–	–
2	Cnd outer1	white/blue	Anode	red
3	Cnd inner2	blue	–	–
4	Cnd outer2/ Shield	black	Shield/GND	green/yellow
5	–	–	Cathode	transparent
6	RTD ref/GND	bare shield	GND/NTC	white & gray
7	RTD sense	red	–	–
8	RTD	green	NTC	green
9	+5V	–	+5V	–

Terminal 4 and 6 are internally connected, either terminal can be used to connect a wire.

* Transparent not connected

TB3 – Conventional (analog) Sensors

Term.	pH		Redox (ORP)	
	Function	Color*	Function	Color
1	Glass	transparent	Platinum	transparent***
2	–	–	–	–
3	Reference**	red	Reference	Shield
4	Solution GND/Shield**	green/yellow and blue	–	–
5	–	–	–	–
6	RTD ref/GND	white	–	–
7	RTD sense	–	–	–
8	RTD	green	–	–
9	+5V	–	–	–

Terminal 4 and 6 are internally connected, either terminal can be used to connect a wire.

* Grey wire not used.

** For pH without solution ground, install jumper 3 to 4.

*** Remove black coating.

TB3 – Conventional (analog) Sensors

Term.	Cond Ind		
	Color. InPro 7250 ST/PFA	Color. InPro 7250 HT	Function
1	Coax inner/transparent	Coax inner/transparent	receive hi
2	red	yellow	receive lo
3	green/yellow	green/yellow	shield/GND
4	brown, white	violet	send lo
5	blue	black	send hi
6	white	white	RTD ref/GND
7	grey	grey	RTD sense
8	green	green	RTD
9	–	–	not used

Terminal 4 and 6 are internally connected, either terminal can be used to connect a wire.

TB3 – Conventional (analog) Sensors

Term.	Dissolved carbon dioxide	
	Color*	Function
1	Coax inner/transparent	Glass
2	–	not used
3**	Coax shield/red	Reference
4**	green/yellow	GND/Shield
5	–	not used
6	white	RTD ref/GND
7	–	not used
8	green	RTD
9	–	+5V

Terminal 4 and 6 are internally connected, either terminal can be used to connect a wire.

* Grey wire not used.

** Jumper 3 to 4 has to be installed.

TB3 – Conventional (analog) Sensors


Term.	Color	Function
1	Coax inner/pink	FET
2	–	not used
3*	yellow	Reference
4*	green / yellow	GND / Shield
5	–	not used
6	white	RTD ref / GND
7	–	not used
8	grey	RTD
9	brown	+5V

Terminal 4 and 6 are internally connected, either terminal can be used to connect a wire.

* Jumper 3 to 4 has to be installed.

4 Wiring example for pH Transmitter (using TB3)

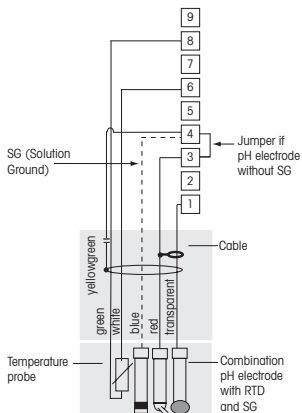pH measurement with monitoring of glass electrode

- 9 +5V
- 8 RTD
- 7 RTD sense
- 6 RTD ref/GND
- 5 not used
- 4 Solution GND/Shield
- 3 Reference
- 2 not used
- 1 Glass

Note: Wire colors only valid for connection with VP cable, grey not connected.

5 General Setup (applies for all parameters) (PATH: Menu/Quick Setup)

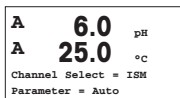
While in Measurement mode press the [MENU] key to bring up the Menu selection. Select "Quick Setup" and press the [ENTER] key.

Display Convention:

1st line on display → a 3rd line on display → c
2nd line on display → b 4th line on display → d

Select the units of measurement for a and b. Only lines a and b can be configured in Quick setup. Go to the Configuration Menu to configure lines c and d.

Channel Selection

Please select the type of Sensor:

Analog: For conventional analog sensors (will be displayed on channel "A").

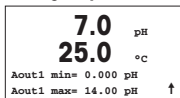
ISM: For ISM sensors (will be displayed on channel "B").

Please select the parameter requirement:

The choice of parameter depends on the level of transmitter. If an ISM sensor is selected, the setting "Auto" means, all possible ISM sensors will be recognized and accepted. If a special parameter is chosen, only this parameter will be recognized and accepted on the transmitter.

**By using analogue 2-e or 4-e conductivity sensors,
see "Conductivity Quick Setup" below for intermediate steps.**

Analog outputs


By selecting Yes the linear 4–20 mA analog output Aout1 will be set up for measurement a when [ENTER] is pressed. Selecting No means that no analog output is set up.

Aout1 min, Aout1 max are the minimum and maximum measurement values for the 4 and 20 mA values respectively.


Note for multi channel devices: The user can configure output A out 3 to measurement c by going back to the previous menu and selecting c.

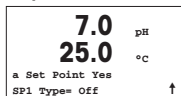Setpoints

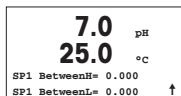
After configuring the Analog Output a setpoint can be configured for that output. If No is selected and [ENTER] is pressed then the quick setup is done and the menus are exited without setting up any setpoint.

Selecting Yes and pressing [ENTER] means a setpoint can be configured for channel a.

Following setpoint Types can be selected:

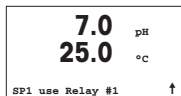- High (High value has to be set)
- Low (Low value has to be set)
- Between (High and Low value has to be set)
- Outside (High and Low value has to be set)

After setting the setpoint value(s) a Relay (none, 1, 2, 3, 4) can be configured for that setpoint. The Relay delay is set to 10 seconds and the Hysteresis is set to 5%. If a relay is configured as CLEAN relay, it will not be selectable in this menu.

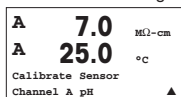
For multi-channel devices: Also setpoints 5 and 6 can be configured and the user can configure a setpoint to measurement c by going back through previous menus and selecting c.

6 pH Calibration incl. ISFET (PATH: Cal)

For Display setup, analog outputs and setpoints see the section “General Setup”.

pH (resp. ISFET) Two point Calibration


Remark: When using ISFET sensors, you have to adjust the zero point first.


While in Measurement mode press the ► key. If the display prompts you to enter the calibration security code, press the ▲ or ▼ key to set the calibration security code, then press the [ENTER] key to confirm the calibration security code.

Press the ▲ or ▼ key to select the pH (resp. ISFET) calibration sub function.

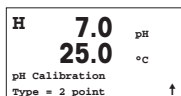
A flashing “H” shows the ongoing calibration process.

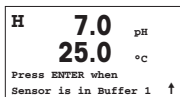Select 2 point Calibration by pressing the [ENTER] key.

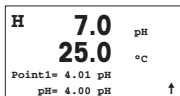
Place the electrode in the first buffer solution and then press the [ENTER] key.


Auto mode: The display shows the buffer that the transmitter has recognized (Point 1) and the measured value.

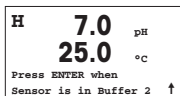Manual mode: Enter the buffer value and press [ENTER] to proceed.

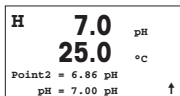
As soon as the stabilisation criteria have been fulfilled (or [ENTER] was pressed in manual mode) the display changes and prompts you to place the electrode in the second buffer solution.


As soon as the stabilisation criteria have been fulfilled (or [ENTER] was pressed in manual mode) the display changes to show the slope calibration factor S and the offset calibration factor Z. Select ADJUST to save the calibration values and the successful Calibration is confirmed on the display. If you select calibrate, the values will not be taken. If an ISM Sensor is used, the Calibration data will be stored in the calibration history.

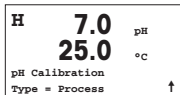Process calibration

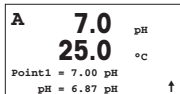Select Process Calibration by pressing the ▲ key once followed by the [ENTER] key. To show the ongoing Calibration Process an "H" is displayed in the top left hand corner.

The "H" changes to "A" (or "B", depends on type of sensor) if Process Calibration is selected to show the user the ongoing calibration on Channel "A" or "B".

Take a sample and press the [ENTER] key again to store the current measuring Value.

After determining the pH Value of the Sample press the ► key again to proceed with the calibration. If the display prompts you to enter the calibration security code, press the ▲ or ▼ key to set the calibration security code, then press the ENTER key to confirm the calibration security code.

A 7.0 pH
25.0 °C
pH S=100.0 % Z=7.124
Save Calibration Yes ↑

Enter the pH value of the sample then press the [ENTER] key to start calibration.

After the calibration the slope calibration factor S and the offset calibration factor Z are displayed. Select ADJUST to save the new calibration values and the successful Calibration is confirmed on the display. If you select calibrate, the values will not be taken. If an ISM Sensor is used, the Calibration data will be stored in the calibration history. The "A" in the top left hand corner disappears.

7 O₂ Calibration (PATH: Cal)

For Display setup, analog outputs and setpoints see the section "General Setup".

O₂ Calibration

A 21.7 pH
A 25.0 °C
Calibrate Sensor
Channel B Oxygen ▲

For multi-channel devices: Using the ▲ and ▼ key on the "Channel A" field lets the user change the channel to be calibrated.

Enter Calibration mode by pressing the ► key.

A flashing "H" (alternating with "A" or "B" to show the channel being calibrated for multi-channel devices) in the top left hand corner shows the ongoing calibration process.

H 21.7 %sat
25.0 °C
Point1 = 100.0 %sat
DO = 0.033 %sat ↑

A DO sensor calibration is always a one point calibration either in Air (Slope) or a zero (Offset) calibration.

It is possible to select a Slope or Zero calibration. A one point slope calibration is done in air and a one point offset calibration is done at 0 ppb DO. Press the [ENTER] key after selecting Slope or Offset.

H 21.7 %sat
25.0 °C
O2 S=0.019nA Z=0.000nA
Save Calibration Yes ↑

Enter the value for Point 1 including a decimal point. DO is the value being measured by the transmitter and sensor in the units set by the user. Press [ENTER] when this value is stable to perform the calibration.

After the calibration the slope calibration factor S and the offset calibration factor Z are displayed.

H 21.7 %sat
A 25.0 °C
Calibration Successful

Select ADJUST to save the calibration values and the successful Calibration is confirmed on the display. If you select calibrate, the values will not be taken. If an ISM Sensor is used, the Calibration data will be stored in the calibration history.

8 CO₂ Calibration (PATH: Cal)

For Display setup, analog outputs and setpoints see the section "General Setup".

CO₂ One point Calibration

A 180.4 hPa
A 26.1 °C
Calibrate Sensor
Channel A CO2

While in measurement mode press the key [CAL]. Press the ▲ or ▼ key to select the Conductivity calibration and press the [ENTER] key to call the sub function.

H 180.4 hPa
A 26.1 °C
CO2 Calibration
Type = 1 Point

Select 1 point and press [ENTER]

H 137.5 hPa
A 26.1 °C
Press ENTER when
Sensor is in Buffer 1

Place the sensor in the buffer solution and press the [ENTER] key to start the calibration.

A 154.5 hPa
A 26.1 °C
A Point1 = 7.00 pH
A CO2 = 7.07 pH

The display shows the buffer the transmitter has recognized (Point 1) and the measured value.

H 154.5 hPa
A 26.1 °C
pH S=100.0% Z=7.048pH
Save Adjust

As soon as the drift conditions have stabilized the display changes to show the slope calibration factor S and the offset calibration factor Z.

After a successful calibration, the calibration values are taken over (Adjust) or were aborted (Calibrate or Abort).

CO₂ Two point Calibration

Start as in chapter CO₂ One point Calibration described.

Select 2 Point calibration.

A **154.5** hPa
A **26.1** °C
CO₂ Calibration
Type = 2 Point ↑

Place the sensor in the first buffer solution and press the [ENTER] key to start the calibration.

H **137.5** hPa
A **26.1** °C
Press ENTER when
Sensor is in Buffer 1 ↑

The display shows the buffer the transmitter has recognized (Point 1) and the measured value.

154.5 hPa
A **26.1** °C
A Point1 = 7.00 pH
A CO₂ = 7.07 pH ↑

As soon as the drift conditions have stabilized, the display changes and prompts you to place the electrode in the second buffer.

A **122.4** hPa
A **26.1** °C
Press ENTER when
Sensor is in Buffer 2 ↑

Place the sensor in the second buffer solution and press the [ENTER] key to go on with the calibration.

2.8 hPa
A **26.1** °C
A Point2 = 9.21 pH ...
A CO₂ = 8.80 pH ↑

The display shows the second buffer the transmitter has recognized (Point 2) and the measured value.

2.8 hPa
A **26.1** °C
pH S=74.21% Z=6.948pH
Save Adjust ↑

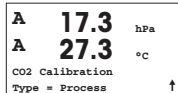
As soon as the drift conditions have stabilized the display changes to show the slope calibration factor S and the offset calibration factor Z.

After a successful calibration, the calibration values are taken over (Adjust) or were aborted (Calibrate or Abort).


CO₂ Process Calibration

Start as in chapter CO₂ One point Calibration described.


Select Process calibration.

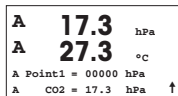
Take a sample and press the [ENTER] key again to store the current measuring value. To show the ongoing calibration process, A or B (depending on the channel) is blinking in the display. After determining the CO₂ value of the sample, press the ► key again to proceed with the calibration.

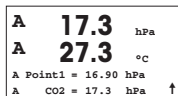Enter the CO₂ value of the sample then press the [ENTER] key to start calibration.

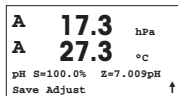The display shows the slope calibration factor S and the offset calibration factor Z. After a successful calibration, the calibration values are taken over (Adjust) or were aborted (Calibrate or Abort).


9 M400 Type 1 Cond Ind only: Conductivity Calibration for inductive Sensors


(PATH: Cal)

For Display setup, analog outputs and setpoints see the section "General Setup".

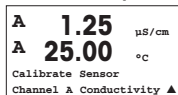
Cond Ind Zero point Calibration

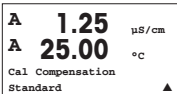
While in measurement mode press the key [CAL]. Press the ▲ or ▼ key to select the Conductivity calibration and press the [ENTER] key to call the sub function.

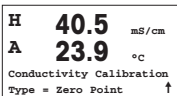Select the kind of calibration (standard, linear or n. water).

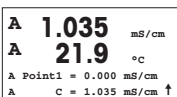Select Zero point and press [ENTER]. Remove the sensor and dry it.


Press the [ENTER] key again to go on with the calibration.

As soon as the conditions have stabilized (or [ENTER] pressed in the manual mode) the display changes to show the multiplier and adder.


After a successful calibration, the calibration values are taken over and stored in the cal history (Adjust), only stored in the cal history (Calibrate) or the calibration can be aborted (Abort).

Cond Ind One point Slope Calibration

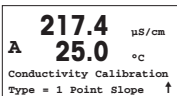Start as in chapter Cond Ind Zero point Calibration described. Select the kind of calibration.

Select 1 point Slope and press [ENTER].

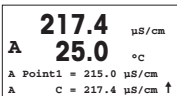Enter the value of calibration point 1 and then press the [ENTER] key to start calibration.

The cell multiplier and the adder are displayed. After a successful calibration, the calibration values are taken over and can be stored in the cal history (Adjust), only stored in the cal history (Calibrate) or the calibration can be aborted (Abort).

10 Conductivity Quick Setup (PATH: Menu/Quick Setup)

For Display setup, analog outputs and setpoints see the section "General Setup".

Sensor Type Selection

1.25 $\mu\text{S}/\text{cm}$
25.0 $^{\circ}\text{C}$
Sensor Type = Cond(2) ▲

Select the type of sensor to be used with the M400 transmitter. Choices are "Cond(2)", used for all 2-Electrode type sensors and "Cond (4)" for all 4-electrode sensors. Press the [ENTER] key.

Cell Constant

1.25 $\mu\text{S}/\text{cm}$
25.0 $^{\circ}\text{C}$
p M=0.1003 A=0.0000
s M=1.0000 A=0.0000

Enter the appropriate cell constant(s):from the sensor label or certificate (M) for 2-electrode sensors, leaving (A) at 0.000; or (M) and (A) values for 4-electrode sensors. Press the [ENTER] key.

Measurement units

1.25 $\mu\text{S}/\text{cm}$
25.0 $^{\circ}\text{C}$
a S/cm
Analog Output? Yes ▲

Select the measurement (conductivity or temperature) and units for measurement. If using analog output, select Yes. Refer back to section "General Setup" to continue setup.

11 Conductivity Calibration for 2-e and 4-e Sensors (PATH: Cal)

For Display setup, analog outputs and setpoints see the section "General Setup".


Cond One point Calibration

A	1.25 $\mu\text{S}/\text{cm}$
A	25.0 $^{\circ}\text{C}$
Calibrate Sensor	
Channel A Conductivity ▲	

While in Measurement mode press the [CAL] key. If the display prompts you to enter the calibration security code, press the ▲ or ▼ key to set the calibration security code, then press the [ENTER] key to confirm the calibration security code.


Press [ENTER] to select the Conductivity calibration sub function.

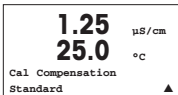A flashing "H" (alternating with "A" or "B" to show the channel being calibrated for multi-channel devices) in the top left hand corner shows the ongoing calibration process.

Choose the compensation mode by using the ▲ or ▼ key.

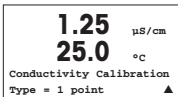
Press [ENTER] to confirm the compensation mode. .

Select "1 point" instead of "2 point Calibration" by pressing the ▼ key once followed by the [ENTER] key.

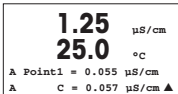
CAUTION: Rinse sensors with a high-purity water solution before every calibration to prevent contamination of the reference solutions.

Place the electrode in the reference solution.

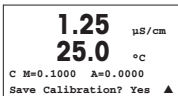Enter the value of Point 1 and press the [ENTER] key.

After the calibration of the cell multiplier or slope calibration factor "M" i.e. cell constant and the Adder or offset calibration factor "A" are displayed.

Select "Yes" to save the calibration values and the successful Calibration is confirmed on the display.

The user gets the message "Re-install sensor" and "Press [ENTER]" on the display. After pressing [ENTER] the M400 returns to the measuring mode.

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© Mettler-Toledo AG 02/2010.
- CH** Mettler-Toledo (Schweiz) GmbH, Im Langacher, Postfach, CH - 8606 Greifensee
- D** Mettler-Toledo GmbH, Prozeßanalytik, Ockerweg 3, D - 35396 Gießen
- F** Mettler-Toledo, Analyse Industrielle S.A.S., 30, Boulevard Douaumont, BP 949, F - 75829 Paris Cedex 17
- CN** Mettler-Toledo Instruments (Shanghai) Co. Ltd., 589 Gui Ping Road, Cao He Jing, CN - 200233 Shanghai
- UK** Mettler-Toledo LTD, 64 Boston Road, Beaumont Leys, GB - Leicester LE4 1AW