

Short instructions AutoRep E™

1. Starting position:
locking lever in upper position
2. To switch on:
press ENTER key or insert
ENCODE Tip and push the locking
lever down to lowest position
3. To adjust volume:
see menu instructions
default mode: DISP (Dispensing)
4. Press STEP key for aspirating and
dispensing liquid
5. Remaining volume in ENCODE Tip
(residual stroke):
discharge by holding down the
STEP key

Operating elements:



- 1 Display
- 2 $\nabla\Delta$ keys (+/-)
- 3 ENTER key
- 4 Menu selection key
- 5 Battery compartment lid
- 6 STEP key
- 7 Locking lever

Menu Instructions

- press MENU key briefly
- hold down MENU key (> 3 s)


Adjusting volume:

MENU 1x \blacktriangleright :



AUTO-DISP-Mode:


MENU 2x \blacktriangleright : $\nabla\Delta$ (ON/OFF)

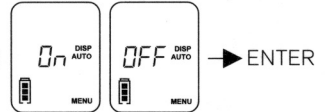

DISP-/PIP-Mode:


MENU 1x $\mathbb{I}\mathbb{I}\mathbb{I}\mathbb{I}\blacktriangleright$ + 1x \blacktriangleright : $\nabla\Delta$ (DISP/PIP)

More information, see operating
manual, pages 22-26.

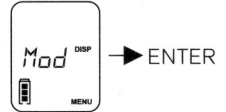


Adjusting speed:

MENU 1x $\mathbb{I}\mathbb{I}\mathbb{I}\mathbb{I}\blacktriangleright$: $\nabla\Delta$ (Aspiration speed)

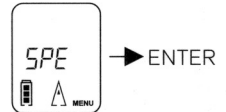

$\nabla\Delta$ (dispensing speed)

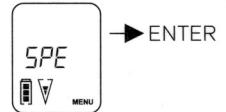

Notes:




Reference run \blacktriangleright press STEP key

Reset:

- remove battery pack or
- simultaneously press the ENTER key and the MENU key.

Additional functions, see operating manual.

Display:

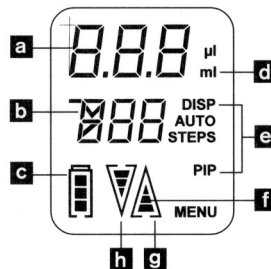

- Tip capacity/Step volume
- Tip/Number of steps
- Battery status
- Volume unit
- Function
- Speed
- Aspirating
- Dispensing