# METTLER TOLEDO DM143-SC

**Double platinum pin electrode**  
**Doppelplatinstift-Elektrode**  
**Electrode à double pointe de platine**  
**Electrodo con doble pin de platino**  
**Двойной платиновый электрод**  
**双铂针电极**

**Double platinum pin electrode**  
for voltametrically and amperometrically  
indicated titrations

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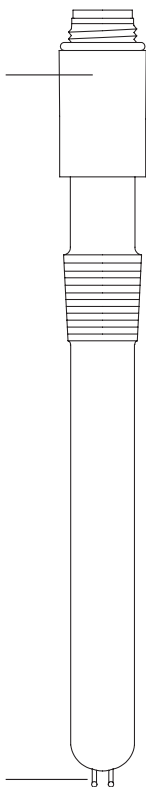
## **1. Pre-operation treatment**

- 1.1 No pre-operation treatment is necessary. The electrode is immediately ready for use.

## **2. Testing**

Only execute testing with a clean and dry electrode, refer to points 4 and 5.

- 2.1 With both electrode poles the resistance between platinum pin and triaxial connector may not exceed  $30 \Omega$  (= electrical continuity).
- 2.2 The resistance between both platinum pins must be greater than  $1 \text{ M} \Omega$  (= insulation).
- 2.3 If the values according to 2.1 and 2.2 cannot be attained, you should replace the electrode.

## **3. Storage and lifetime**

- 3.1 The electrode should be stored clean and dry.
- 3.2 The lifetime of the electrode, depending on its proper use and treatment, is almost unrestricted.

## **4. Maintenance Instructions**

- 4.1 Polarizing currents which are too high cause an electrochemical contamination of the platinum pins.
- 4.2 Do not let the triaxial connector corrode.

## **5. Cleaning procedures**

- 5.1 Place the electrode in an ultrasonic bath with deion. water or ethanol for a few minutes.
- 5.2 Place the electrode in sulfochromic acid for 60 s and rinse it thoroughly with deion. water and ethanol.

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