METTLER TOLEDO DM140-SC

Platinum ring electrode
Platinring-Elektrode
Electrode à l'anneau de platine
Electrodo con anillo de platino
Электрод с платиновым кольцом
金属铂环电极

Combined platinum ring electrode
with ceramic frit for redox titrations.

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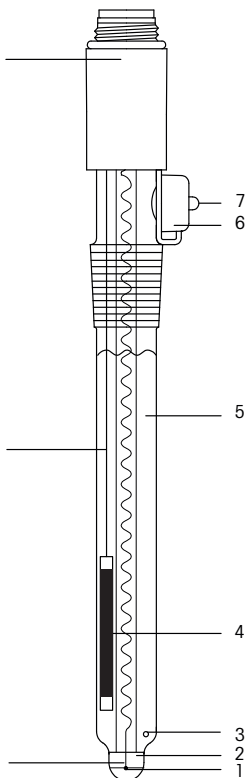
1. Start-up maintenance

- 1.1 To equalize pressure remove the rubber cap (6) or perforate its nipple (7).
- 1.2 Refill the reference electrolyte (5) – (e.g. METTLER TOLEDO No. 51340049) – to approx. 1 cm below the aperture (6).
- 1.3 Place the electrode in deion. water for 15 minutes to dissolve the KCl crystals in or on the frit (3).

2. Testing

The electrode can be tested with a redox buffer solution, i.e. its potential and response time are measured.

- 2.1 The electrode potential should correspond to the values indicated on the buffer solution (e.g. METTLER TOLEDO No. 51340065 or 5134006).
- 2.2 Response time of the electrode: After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 2.3 Testing a used electrode, a comparison between titration curves is suitable: perform a titration with sodium thiosulfate (0.1 mol/L) and iodine (0.1 mol/L) with the new electrode. The titration goes fast with a big steep potential change at the equivalence point. A used electrode performing the same titration should present both these parameters.

3. Storage and lifetime

- 3.1 Store the electrode in the reference electrolyte – the frit has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 3.2 Never let the electrode go dry, otherwise the KCl crystals in or on the frit have to be dissolved.
- 3.3 The lifetime of the electrode is, depending on its proper use and treatment, almost unrestricted.

4. Maintenance instructions

- 4.1 The tolerable temperature range extends from 0 to 80 °C (32 – 176 °F).
- 4.2 In strong oxidizing solutions an oxygen adsorption takes place on the platinum surface of the electrode, in strong reducing solutions hydrogen adsorption. This can cause a change in the redox potential as well as an increased response time.
- 4.3 The sample solution should never enter into the reference electrode through the frit. Therefore the level of the reference electrolyte must be always higher than the sample solution level.

- 4.4 No reaction should take place between the sample solution and the reference electrolyte: sulfides, bromides, iodides and cyanides react with Ag^+ and some cations (Ag^+ , Hg^{2+} , Au^{3+} , Pb^{2+}) react with Cl^- of the reference electrolyte and cause a blockage of the frit.
- 4.5 Solutions containing protein quickly block the frit, due to protein precipitates.
- 4.6 Do not let the cable connector plug corrode.

5. Cleaning procedures

- 5.1 Contamination of the electrolyte: drain the electrolyte solution and refill with new electrolyte. Place the electrode in deion. water for several hours.
- 5.2 Frit blocked by AgCl : place it overnight in concentrated ammonia, rinse and condition it for 1 hour in a buffer solution pH 4.
- 5.3 Frit blocked by sulfide precipitate: place it in a solution of thiourea (8% thiourea /1 mol/ L HCl, e. g. METTLER TOLEDO No. 51340070) until the dark color has vanished.
- 5.4 Frit blocked by protein: place it for at least 1 hour in pepsin solution (5% pepsin /0.1 mol/L HCl, e.g. METTLER TOLEDO No. 51340068).
- 5.4 Frit blocked by protein: place it for at least 1 hour in pepsin solution (5% pepsin/0.1 mol/L HCl, e.g. METTLER TOLEDO No. 51340068).
- 5.5 Contamination of the platinum surface: depending on the contamination several cleaning procedures can be carried out.
 - a) Place the electrode for a few minutes in an ultrasonic bath with deion. water or ethanol.
 - b) Place the electrode for one minute in chromium sulfuric acid or aqua regia and rinse it thoroughly with deion. water.
 - c) Place the electrode for 30 minutes in a pH 4 buffer solution (Potassium hydrogen phthalate) saturated with quinhydrone to remove the absorbed oxygen or hydrogen.
 - d) Polish the platinum surface with Al_2O_3 powder.

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