Quick Setup Guide **InPro 8300 RAMS**

For more detailed information please consult the instruction manual.

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METTLER TOLEDO

1 Contents of package

The package contains the following components:

- InPro 8300 RAMS
- Screwdriver
- Face spanner
- Installation wrench for OPL bits
- 2 OPL bits
- Container with spare desiccant capsules
- 2 × O-rings, 34.59 × 2.62
- Manual
- Software CD (for types InPro 8300 RAMS COMBINE and CALI only)

2 Installation of the InPro 8300 RAMS

2.1 Mechanical installation

The RAMS can be fitted to existing VARINLINE® measuring cells or VARINLINE® sight glasses. Make sure that the measuring cells are welded into the pipe so that the electronic parts of the InPro 8300 RAMS face each other horizontally.

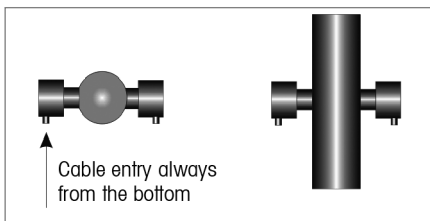
Fig. 1: Installation position of the measuring cell

Insert the OPL bits in the VARINLINE® adapters in the measuring cell and secure them using hinged clamps; tool required: a/f 10 mm open-ended, ring or socket spanner. Make sure that the recesses for the securing pins (Fig. 2, item 16) are facing downwards. This way you ensure that the cable glands will also point downwards once you have fitted the RAMS to the pipe.

Carry out the following steps for both electronic parts:

- Remove the protective caps from the InPro 8300 RAMS OPL-bits (Fig. 2, item 7).
- Insert a 34.59 × 2.62 mm O-ring into the groove (Fig. 2, 18).
- Take the desiccant capsule out of the tube and place it in the recess (Fig. 2, item 17) in the electronic part. Immediately close the desiccant tube again!
- Fit the electronic part to the VARINLINE® adapter or the OPL bit and use the face spanner to fasten it. The active electronic part (Fig. 2, item 4) has to be mounted at the side with the shorter OPL bit.

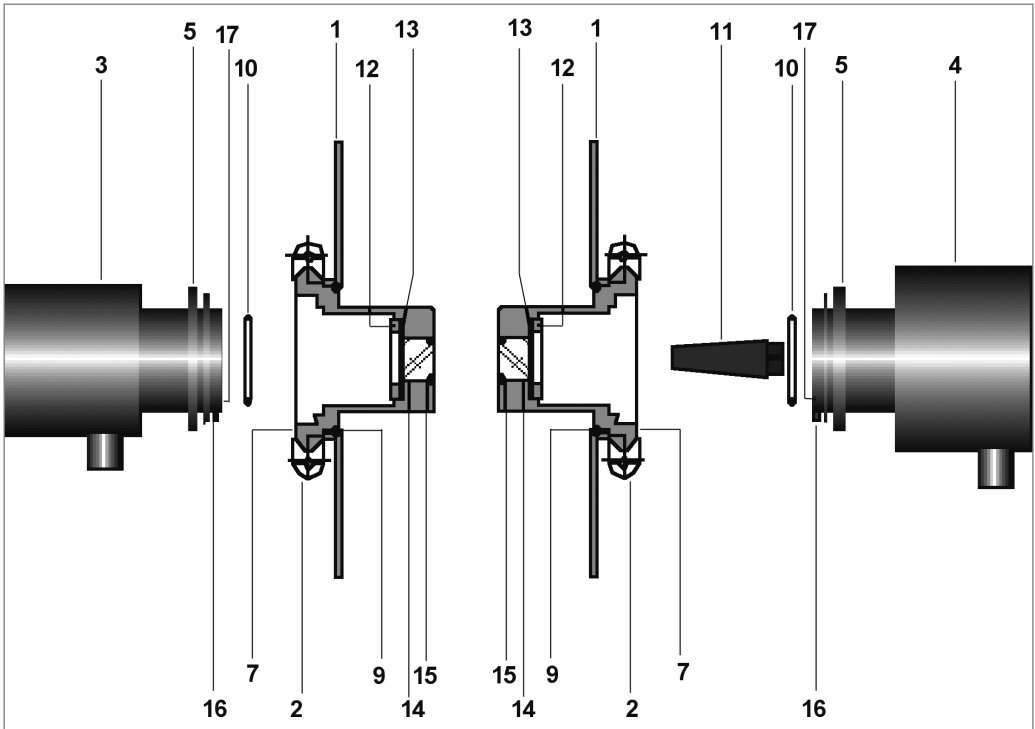
Fig. 2

- 1 VARINLINE® housing (in pipeline)
- 2 Hinged clamps
- 3 Passive electronic part of the InPro 8300 RAMS
- 4 Active electronic part of the InPro 8300 RAMS (when using an OPL bit and a 68 mm window always mount it on the side where the big window is fitted, when using 2 OPL bits mount it on the side where the short OPL bit is fitted)
- 5 Threaded ring
- 7 OPL bit
- 9 O-ring, 60 × 3
- 10 O-ring, 34.59 × 2.62
- 11 Spacer (only when 2 OPL bits are used)
- 12 Threaded ring
- 13 Thrust ring
- 14 Window, 28 mm
- 15 O-ring, 22 × 3
- 16 Securing pin
- 17 Recess for desiccant capsule

2.1 Electrical installation

The InPro 8300 RAMS is supplied with the connection cable firmly connected between active and passive part. Once the InPro 8300 RAMS has been installed mechanically, the electrical connections can be made.

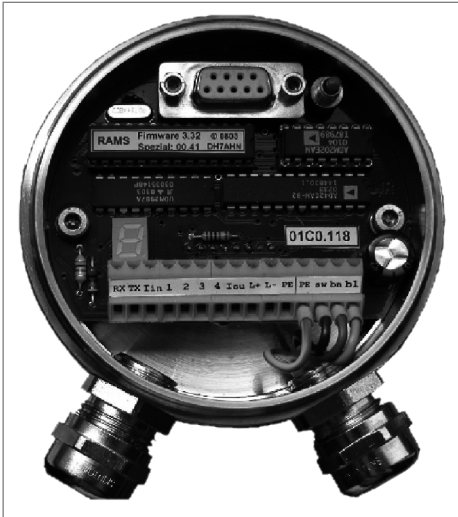
Fig. 3

- Unscrew the window plate from the InPro 8300 RAMS
- Pass the stripped cable through the cable gland
Cable diameter 4 ... 10 mm,
Core cross section 0.75 mm²
- Connect the cable cores in accordance with the circuit diagram
- ⚠ **IMPORTANT:** Input voltage is 24 VDC, 200 mA only
- Insert the desiccant capsule in the terminal compartment
- Close the window plate

Tools required:

- Open-ended wrench d/f 22 mm (not included)
- Cable stripper (not included)
- Screwdriver 2.4 mm (included in the InPro 8300 RAMS accessories kit)
- Desiccant capsules (2 pcs, included in the InPro 8300 RAMS accessories kit)

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- CH** Mettler-Toledo (Schweiz) GmbH, Im Langacher, Postfach, CH - 8606 Greifensee
- D** Mettler-Toledo GmbH, Prozeßanalytik, Ockerweg 3, D - 35396 Gießen
- F** Mettler-Toledo, Analyse Industrielle S.A.S., 30, Boulevard Douaumont, BP 949,
- CN** Mettler-Toledo Instruments (Shanghai) Co. Ltd., 589 Gui Ping Road, Cao He Jing,
- UK** Mettler-Toledo LTD, 64 Boston Road, Beaumont Leys, GB - Leicester LE4 1AW
- USA** Mettler-Toledo, Process Analytics, Inc., 36 Middlesex Turnpike, Bedford, MA 01730, USA