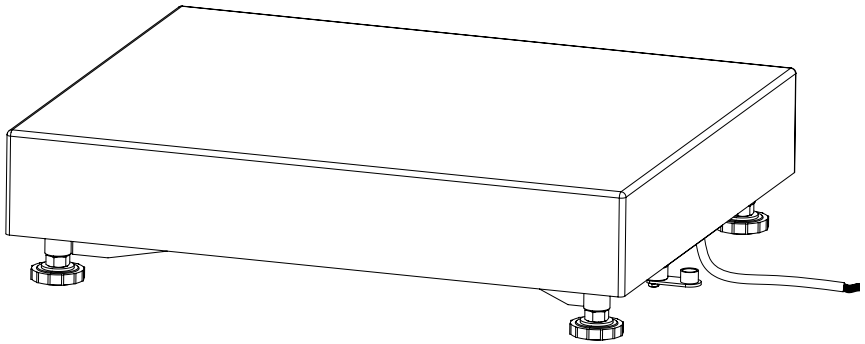
# PBA426 / PBA426x

## Weighing Platform

**Bedienungsanleitung** Wägeplattform **PBA426 / PBA426x**  
**Manuel Utilisateur** Plate-forme de pesage **PBA426 / PBA426x**  
**Manual de usuario** Plataforma de pesaje **PBA426 / PBA426x**  
**Manuale utente** Piattaforma di pesata **PBA426 / PBA426x**  
**Manual do usuário** Plataforma de pesagem **PBA426 / PBA426x**

METTLER TOLEDO

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## ServiceXXL

Tailored Services

Congratulations on choosing the quality and precision of METTLER TOLEDO. Proper use according to these instructions and regular calibration and maintenance by our factory-trained service team ensure dependable and accurate operation, protecting your investment. Contact us about a ServiceXXL agreement tailored to your needs and budget.

We invite you to register your product at

[www.mt.com/productregistration](http://www.mt.com/productregistration)

so we can contact you about enhancements, updates and important notifications concerning your METTLER TOLEDO product.

# Contents

<b>1</b>	<b>Safety Instructions</b> .....	<b>6</b>
<b>2</b>	<b>Setting up the Weighing Platform</b> .....	<b>7</b>
2.1	Selecting the site of installation.....	7
2.2	Leveling .....	8
2.3	Connecting the weighing platform to the weighing terminal .....	9
2.4	Equipotential bonding (PBA426x).....	9
<b>3</b>	<b>Operation Limits</b> .....	<b>10</b>
<b>4</b>	<b>Cleaning</b> .....	<b>11</b>
<b>5</b>	<b>Technical Data</b> .....	<b>12</b>
5.1	Technical data of the weighing cell.....	12
5.2	Dimensions.....	12
5.3	Accessories .....	13
<b>6</b>	<b>Appendix</b> .....	<b>14</b>
6.1	Notice for verified weighing instruments in EC countries .....	14
6.2	Disposal .....	14
6.3	Declaration of conformity .....	15

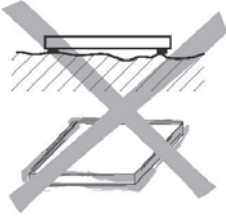
# 1 Safety Instructions

Product safety plays an important role at METTLER TOLEDO. Non-observance of the following instructions can lead to damage to the weighing platform and/or injuries.

- ▲ Read this manual carefully before operating or servicing this equipment.
- ▲ Strictly observe these instructions and save this manual for future reference.
- ▲ Only personnel trained and qualified by METTLER TOLEDO may install and maintain the weighing platform.
- ▲ Be careful when transporting or lifting heavy devices.
- ▲ Always disconnect this device from the power source before installing, servicing, cleaning or performing maintenance.
- ▲ Make sure the weighing platform reached room temperature before switching on the power supply.
- ▲ The explosion-protected weighing platforms (PBA426x) are approved for operation in following hazardous areas: ATEX Category 2, gas/dust (Zone 1/21) - FM I.S. Class I, II, III Div. 1 (Load Cell FM approved)
- ▲ There is an increased risk of injury and damage when the weighing platforms are used in hazardous areas! Special care must be taken when working in such areas.
- ▲ Explosion-protected weighing platforms may only be used in hazardous areas in conjunction with weighing terminals with the appropriate approval and interface specification.
- ▲ The connection cable may not be disconnected from the weighing terminal while energized.
- ▲ In case of PBA426x only use cable glands that are suitable and approved for hazardous areas for inserting the weighing cell cable into the service terminal.

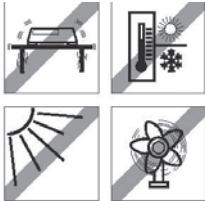
## 2 Setting up the Weighing Platform

### 2.1

#### Selecting the Site of Installation

- The surface must be able to bear the weighing platform under maximum load at the points of support. At the same time it should be so stable that no vibrations arise during weighing. This is also to be observed when installing the weighing platform in conveyor and similar systems.
- If possible, platform installation should be located in an area free from vibration.

#### Ambient Conditions

- Observe the following ambient conditions:
  - No direct sunshine
  - No permanent water on the load cell (max. 120 min. per day)
  - Absolute humidity shall stay below 30 g/m<sup>3</sup> (environment)
  - No excessive temperature fluctuations
  - Temperature range –10 °C to +40 °C

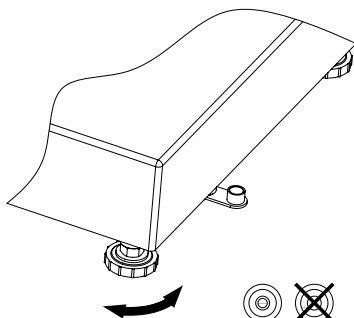
## 2.2

### Leveling

Only a weighing platform which is aligned exactly horizontally supplies exact weighing results.

The weighing platform has to be leveled during the initial installation and whenever its location is changed.

1. Turn the adjustable feet of the weighing platform until the air bubble of the spirit level is positioned in the inner circle.
2. Tighten the lock nuts of the adjustable feet.

## 2.3 Connecting the Weighing Platform to the Weighing Terminal

The weighing platform PBA426(x) is designed for use with analog METTLER TOLEDO weighing terminals.

1. Insert the weighing platform cable through the cable gland into the weighing terminal.
2. Connect the weighing platform cable to the 7-pin terminal strip of the weighing terminal according to the following tables.

PBA426-... / PBA426x-...	
Terminal	Color
EXC+	blue
EXC-	black
SIG+	white
SIG-	red
SEN+	green
SEN-	grey
Shield	yellow

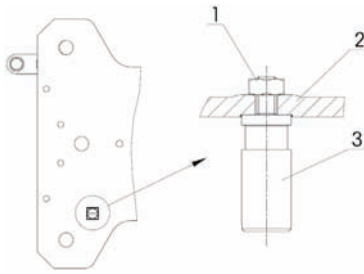
## 2.4 Equipotential Bonding (PBA426x)

The equipotential bonding for the PBA426x platform series is mounted in the factory.

The wiring must be installed by a professional electrician on customer site. METTLER TOLEDO Service only has a monitoring and consulting function here.

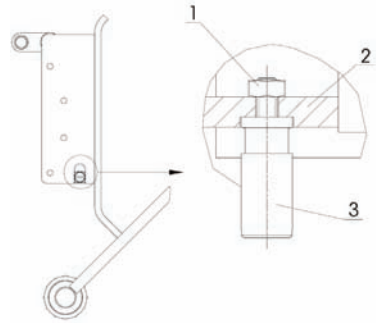
Connect equipotential bonding (PA) of all devices (weighing platform and service terminal) in accordance with the country-specific regulations and standards. In the process, make sure that all device housings are connected to the same potential via the PA terminal.

Note: The equipment is protected against electric shock acc. IEC60950, class II.

**PBA426x - A/QA**

- (1) Hexagonal lock nut M4 KEPS (119319)
- (2) Base frame
- (3) Equipotential bonding terminal (72200560)

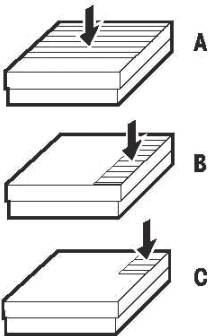
**PBA426x - BB/B/QB**

- (1) Hexagonal lock nut M4 KEPS (119319)
- (2) Base frame
- (3) Equipotential bonding terminal (72200560)

### 3 Operation Limits

With the robust scale design you can occasionally exceed the rated capacity of the platform without damage. The maximum static safe load must never be exceeded.

→ Observe the following operation limits.

Model	A – center load	B – side load	C – corner load
PBA426(x)-A...	40 kg / 80 lb	30 kg / 60 lb	15 kg / 30 lb
PBA426(x)-BB...	100 kg / 200 lb	70 kg / 140 lb	35 kg / 70 lb
PBA426(x)-B...	200 kg / 400 lb	140 kg / 280 lb	75 kg / 150 lb
PBA426(x)-QA...	40 kg / 80 lb	30 kg / 60 lb	15 kg / 30 lb
PBA426(x)-QB...	100 kg / 200 lb	70 kg / 140 lb	35 kg / 70 lb

→ Avoid falling loads, shock loads as well as impacts from the side.

## 4 Cleaning

- Take off the load panel and remove any dirt and foreign substances which may have collected underneath it. Do not use any hard objects to do so. Do not disassemble the weighing platform.
- Only use disinfectants and cleaning agents in accordance with the manufacturers's specifications and instructions and then rinse with clear water.
- Clean the weighing platform with splash water (12.5 L/min.) and remove any remaining detergent by rinsing with clear water.
- To prolonge the life time of the load cell, dry it off with a soft lintfree cloth immediately after cleaning.

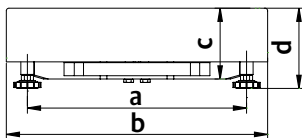
## 5 Technical Data

### 5.1 Technical Data of the Weighing Cell SLP535 (PW15)

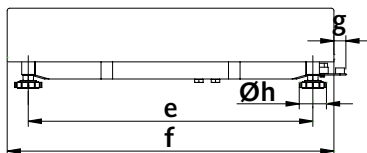
Sensitivity	2 mV/V +/- 0.2 mV/V
Input resistance	380 $\Omega$ +/- 15 $\Omega$
Output resistance	359 $\Omega$ +/- 5 $\Omega$
Supply voltage	Recommended: 5 V DC/AC Maximum: 12 V DC/AC
Approval	OIML / NTEP / ATEX / FM
Ignition protection type	II2G Ex ia IIC T4/T6 -30°C<T <sub>a</sub> <40°C(T6)70°C(T4) II2D Ex ia D21 T100°C
IP protection type	IP65/IP67

### 5.2 Dimensions

Front view

Side view

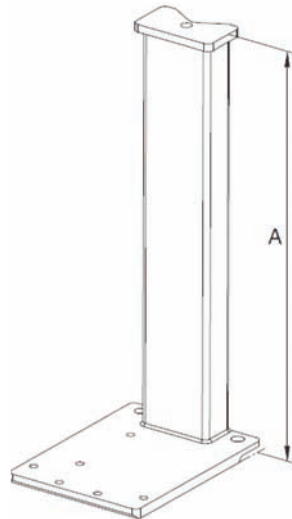
Model – Dimension in mm / "	a	b	c	d*	e	f	g	h
PBA426(x)-A... mm	175	240	59	86	235	300	21	42
" "	6,89	9,45	2,32	3,38	9,25	7,87	0,83	1,65
PBA426(x)-BB... mm	235	300	76	97	335	400	18	42
" "	9,25	11,81	2,99	3,81	13,18	15,74	0,7	1,65
PBA426(x)-B... mm	335	400	108,5	123	435	500	18	42
" "	13,18	15,74	4,27	4,84	17,12	19,68	0,7	1,65
PBA426(x)-QA... mm	163	228	59	86	163	228	21	42
" "	6,41	8,97	2,32	3,38	6,41	8,97	0,83	1,65
PBA426(x)-QB... mm	240	305	76	97	254	305	17	42
" "	9,45	12,0	2,99	3,81	10,0	12,0	0,67	1,65

\*with the hygienic Kit Option, the platform heights (d) increases by 11 mm / 0.43"

### 5.3

### Accessories

Designation	Order No.
Stand (column) A terminal specific column adapter might be needed. The order number can be found in the user manual for the terminal.	
Stainless steel, 120 mm / 4.7" high	72 229 393
Stainless steel, 330 mm / 13" high	72 198 702
Stainless steel, 660 mm / 26" high	72 198 703
Stainless steel, 900 mm / 35.4" high	72 198 704

**For installation instructions please see installation manual no. 72 205 196A**

Note: For PBA426x only these columns might be used in combination with the platform (stainless steel 1.4301)

## 6 Appendix

### 6.1

#### Notice for verified Weighing Instruments in EC Countries

Weighing instruments verified at the place of manufacture bear the preceding mark on the packing label and a green "M" sticker on the descriptive plate. They may be set to work immediately.

Weighing instruments which are verified in two steps have no green "M" on the descriptive plate and bear the preceding identification mark on the packing label. The second step of the verification must be carried

out by the approved METTLER TOLEDO service or by the W & M authorities. Please contact your METTLER TOLEDO organization. The first step of the verification has been carried out at the manufacturing plant. It comprises all tests according to EN45501-8.2.2.

If national regulations in individual countries limit the period of validity of the certification, the operator of such a scale is himself responsible for its timely re-certification.

### 6.2

#### Disposal

In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE), this device must not be disposed of in domestic waste. This also applies to countries outside the EU as per their specific regulations.

→ Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.

If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

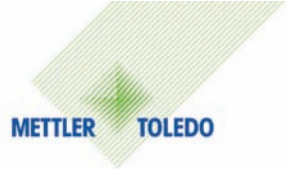
Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related. Thank you for your contribution to environmental protection.

## 6.3 Declaration of Conformity

# METTLER TOLEDO

Legal Metrology

**Declaration of Conformity**  
Konformitätserklärung  
Déclaration de conformité  
Declaración de Conformidad  
Conformiteitsverklaring  
Dichiarazione di conformità

We, Wir, Nous, Nosotros, Wij , Noi

**Mettler-Toledo (ChangZhou) Measurement Technology Ltd.**  
111 West TaiHu Road, XinBei District, ChangZhou, JiangSu, 213125, P.R.China

**Declare under our sole responsibility that the product,**  
erklären, in alleiniger Verantwortung, dass dieses Produkt,  
déclarons sous notre seule responsabilité que le produit,  
declaramos, bajo nuestra sola responsabilidad, que el producto,  
verklaren onder onze verantwoordelijkheid, dat het product,  
dichiariamo sotto nostra unica responsabilità, che il prodotto,

**Model/Type: PBA426(x) series weighing platform**

**to which this declaration relates, is in conformity with the following standard(s) or other normative document(s).**  
auf das sich diese Erklärung bezieht, mit der/den folgenden Norm(en) oder Richtlinie(n) übereinstimmt.  
Auquel se réfère cette déclaration est conforme à la (aux) norme(s) ou au(x) document(s) normatif(s).  
Al que se refiere esta declaración es conforme a la(s) norma(s) u otro(s) documento(s) normativo(s).  
Waarnaar deze verklaring verwijst, aan de volende norm(en) of richtlijn(en) beantwoordt.  
A cui si riferisce questa dichiarazione è conforme alla/e sequente/i norma/e o documento/i normativo/i.

EC marking	EC Directive:	Applicable Standards.
	94/9/EC ATEX (only for PBA426x)	EN 13463-1 EN 60079-0: 2006 EN 60079-11: 2007 EN 61241-0: 2006 EN 61241-11: 2004
For non-automatic weighing instrument used in an <b>Article 1,2.(a)</b> application ,additional metrological marking according to Annex IV of Council Directive 2009/23/EC must be attached to the instrument.		
	2009/23/EC Non-automatic weighing instruments	EN 45501:1992 <sup>1)</sup>

1) **Applies to certified non-automatic weighing instruments only in connection with an approved and compatible weighing indicator.**

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Issued on: 2010-1-29  
Mettler-Toledo (ChangZhou) Measurement Technology Ltd.

Zhu Dan  
General Manager

Yang JiaWu  
QA Manager

**Important notice concerning verified instruments see overleaf!**

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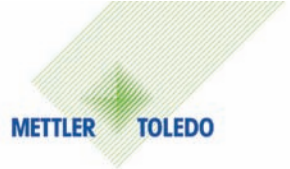
We, Wir, Nous, Nosotros, Wij , Noi

Mettler-Toledo (ChangZhou) Measurement Technology Ltd.  
111 West TaiHu Road, XinBei District, ChangZhou, JiangSu, 213125, P.R.China

Declare under our sole responsibility that the product, erklären, in alleiniger Verantwortung, dass dieses Produkt, déclarons sous notre seule responsabilité que le produit, declaramos, bajo nuestra sola responsabilidad, que el producto, verklaren onder onze verantwoordelijkheid, dat het product, dichiariamo sotto nostra unica responsabilità, che il prodotto,

Model/Type: PBA426(x) series weighing platform

to which this declaration relates, is in conformity with the following standard(s) or other normative document(s). auf das sich diese Erklärung bezieht, mit der/den folgenden Norm(en) oder Richtlinie(n) übereinstimmt. Auquel se réfère cette déclaration est conforme à la (aux) norme(s) ou au(x) document(s) normatif(s). Al que se refiere esta declaración es conforme a la(s) norma(s) u otro(s) documento(s) normativo(s). Waarnaar deze verklaring verwijst, aan de volende norm(en) of richtlijn(en) beantwoordt. A cui si riferisce questa dichiarazione è conforme alla/e sequente/i norma/e o documento/i normativo/i.

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- 1) Applies to certified non-automatic weighing instruments only in connection with an approved and compatible weighing indicator.

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Issued on: 2010-1-29

Mettler-Toledo (ChangZhou) Measurement Technology Ltd.

Zhu Dan  
General Manager

Yang JiaWu  
QA Manager

Important notice concerning verified instruments see overleaf!

[www.mt.com/service](http://www.mt.com/service)

For more information  
Weitere Informationen erhalten Sie unter:  
Pour plus d'information  
Para más información  
Per ulteriori informazioni  
Para mais informações

**Mettler-Toledo AG**  
**CH-8606 Greifensee**  
**Switzerland**  
**Tel. + 41 44 944 22 11**  
**Fax + 41 44 944 30 60**

Subject to technical changes  
© 07/2011 Mettler-Toledo AG  
Printed in Switzerland  
Order number 72230073A

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